

# Nitrogen Fill Movie

This document is recorded as a QuickTime Movie File. It shows how to refill a magnet's liquid nitrogen vessel and gives safety advices.

## **WARNING:**

When this film was recorded some time ago, latex transfer hoses were in use. It has been reported that latex hoses may split during use, therefore it is **IMPERATIVE** that you **do not use latex** as the material for the transfer hose. Hoses with metal sleeving, teflon hoses or surrogated PFA hoses are acceptable for use.

## **SGI systems**

To view the movie on an SGI system, type

```
movieplayer /CDROM/misc/n2fills.mov
```

in a UNIX shell.

## **WINDOWS**

To start the movie under WINDOWS, just click on this line or on the picture below. Install QuickTime from [www.apple.com/quicktime/download/](http://www.apple.com/quicktime/download/) if required.

This movie was prepared by BRUKER INSTRUMENTS INC. Manning Park, Billerica, M.A. All rights are reserved.

Document Part Number: **Z31338**

Drawing Number: **1047001**