How to recover your data

Your data acquired on any of the NMR Service spectrometers are available at all times and from anywhere through our data server. To recover them, you will have to download and install the program [FileZilla](#), and configure it to connect to our server securely.

In order to preserve the privacy of the groups' research data, each research group and company has a virtual account for its exclusive use, that allows it to access only its data folders on the SeRMN spectrometers and computers.

The data server uses [ProFTPd](#) to manage the secure connection and transfer of data using the [W \] protocol FTPS/FTPES](#). This data transfer protocol uses [W TLS](#) to encrypt communications between the client and the server. This is the same encryption protocol that browsers use to access secure web servers via HTTPS.

To access the data server you will need a [W FTP client](#) that allows TLS encrypted connections. We recommend the program [FileZilla](#) although you can use others that support the FTPS/FTPES (FTP over TLS) protocol.

The folders on the data server

When you log in, you will see several folders that link directly to data folders on currently available spectrometers as well as former spectrometers. The name of the folder indicates which spectrometer it corresponds to.

Current spectrometers

These folders correspond to the spectrometers currently installed in the service:

Folder name	Current spectrometers
av-dpx-250-auto	Spectrometer Avance DPX 250 MHz, with sample changer.
av-neo-300	Spectrometer Avance Neo 300 MHz for self-service.
av3-400-sb	Spectrometer Avance III 400 MHz, with an HR-MAS accessory.
av-neo-400	Spectrometer Avance Neo 400 MHz for self-service.
av-neo-500	Spectrometer Avance Neo 500 MHz, with cryoprobe.
av3-600	Spectrometer Avance III+ 600 MHz, with sample changer.
biospec-7T	Spectrometer BioSpec 70/30USR with an horizontal magnet.

Former spectrometers

These folders correspond to spectrometers previously installed in our facility that no longer exist:

Folder name	Former spectrometers
av-dpx-250-robot	Spectrometer Avance DPX-250 MHz, with sample changer.
av-dpx-360	Spectrometer Avance DPX-360 MHz for self-service.

Folder name	Former spectrometers
av2-400-wb	Spectrometer Avancell 400 MHz, wide-bore magnet, with CP/MAS accessory.
av-drx-500	Spectrometer Avance DRX-500 MHz, with cryoprobe and LC-DAD-(SPE)-NMR/MS.

You will probably never see all of these instruments, as each group only sees the spectrometers that at least one member of the group has access to.

Each research group and company only has access to their data folders.

Thus, when acquiring spectra **make sure you save the data in your group folder**. If you save it in the wrong folder, you will not be able to retrieve it through the data server.

Use only unaccented letters, numbers, and dashes (uppercase and lowercase) for experiment names,

It is highly recommended that you only use:

- lowercase and uppercase letters without accents: a-z A-Z
- numbers: 0-9
- hyphens: - _

for experiment names, because other characters (accents, asterisk, quotes, parentheses, etc.) could cause problems when retrieving your data.

Open the firewall of your institute or company

The FTPS/FTPES client and server communicate via ports 990 (FTPS) and 60000-65534 (IANA Ephemeral ports), thus you will have to make sure they are not blocked by any firewalls.

In mid-2022 **we changed the FTPS server port number from 21 to 990** in order to use the official port number for this protocol:

Port 990 Details

Service: ftps

Details: FTPS Protocol, FTP over TLS/SSL (IANA official) uses ports 989 and 990.

<https://www.speedguide.net/port.php?port=990>

If you access our server from within the UAB, or from a research institute or company on campus that has its network integrated into the UAB network, you should have no problems connecting and downloading your data.

Conversely, if you connect from an off-campus computer, and your institute or company has a firewall to protect its network, you will have to ask the people responsible for the firewall to configure it so that you can recover your data .

The configuration is very simple and does not entail an unacceptable security risk, the steps required are :

- open outbound **port 990 (FTPS)** for the **TCP** protocol; and
- open inbound the **range of ports 60000-65534** (IANA Ephemeral ports for FTPS) for the **TCP protocol** to computers on your network


so that you can connect to our FTPS-TLS server via **FTPS-Passive** from inside the firewall.

If they want to further restrict open ports only to our IP address, tell them it's **IP: 158.109.58.175** with **hostname: sermn.uab.cat**

Don't worry if you don't understand this technical information, it is meant for your firewall administrators.

Access your data with FileZilla

The program we recommend to access our data server is [FileZilla](#), an FTP client with a graphical interface, available for Windows, Mac OS X and Linux.

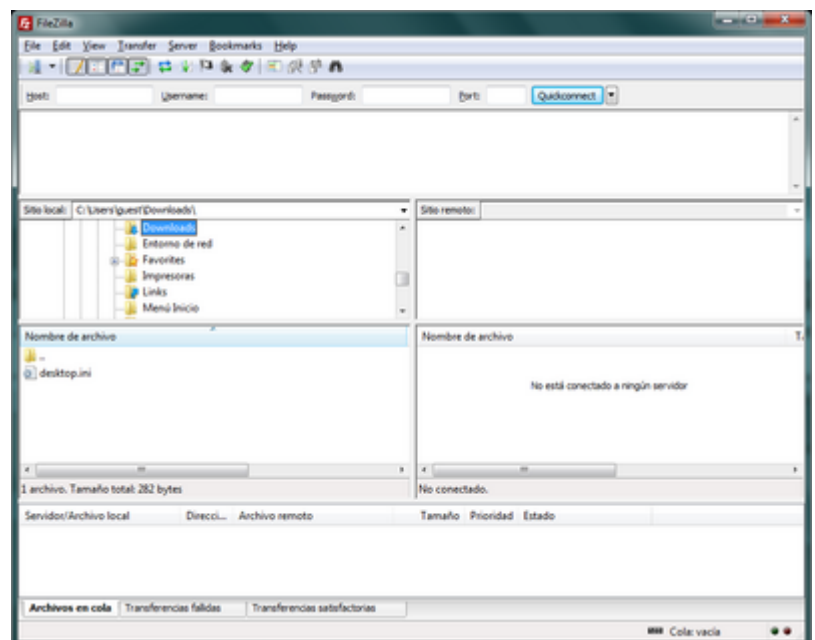At the time of writing these instructions, the most current version is 3.66.4, but since new versions come out quite often, you will probably find a newer version to download and install. On the other hand, it is highly recommended that you pay attention to the program when it informs you that there is a new version and suggests that you download and install it. Doing so is only takes a moment, and requires no software reconfiguration.

Installation

On the website of the FileZilla program you will find [complete, detailed and up-to-date installation instructions](#) on how to install FileZilla Client on MS-Windows, Mac OS X, and GNU/Linux.

You can download the program directly from the [FileZilla website](#).

Detailed configuration

If you are not familiar with FileZilla configuration [follow these detailed instructions](#). There we explain step by step how to configure the program to access our data server, in addition, the multiple screenshots will help you to successfully complete the configuration process.

Otherwise, use the quick setup instructions in the next section.

Quick setup

If you are familiar with FileZilla use these short instructions to set it up:

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