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**2017-05-18 - TREBALL EN MARXA / WORK IN PROGRESS**

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- [Debian multiarch support](#)
- [Multiarch HOWTO for users](#). User-oriented info on how to use multiarch: configuration, installation, cross-building.

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- Install Bruker NMR Software
- Install GPL licensed GNU GCC 4.9.1 compiler (recommended)

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```
Note: Using init directory /etc/init.d  
Note: Using Systemd for service control  
/usr/diskless/prog/bin/setfirewall.sh: install firewall rules for  
Spectrometer Operating System (show only)  
command: iptables-restore < /etc/sysconfig/iptables  
saving of iptables not necessary  
command: ip6tables-restore < /etc/sysconfig/ip6tables  
saving of ip6tables not necessary  
/usr/diskless/prog/bin/setfirewall.sh: terminated with status OK
```

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```
*** Error ***  
  
Could not show firewall rules for MICS:  
(command: /bin/bash /opt/Bruker/mics/setfirewall.sh show)  
  
/opt/Bruker/mics/service-lib.sh: line 241: rpm: command not found  
Fatal error: is not a number
```

```
/opt/Bruker/mics/setfirewall.sh: terminated with status 3
```

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```
*** Warning ***
```

```
Cannot create user account 'flexlm' and/or group 'flexlm':  
Cannot create user account flexlm: useradd flexlm : useradd: group flexlm  
exists - if you want to add this user to that group, use -g.
```

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```
installation of DISKLESS 3.5 finished>  
>> installnmr terminated at 2017-05-18 11:51:47 +0200  
>> STEP post terminated for mod=hostacq with result = NEXT  
>> STEP post started for mod=ipso at 2017-05-18 11:51:47 +0200 ...  
>> cd /tftpboot  
>> Status: Running post-installation scripts for IpsoOS  
>> ::file::write (version.txt, -log)  
>> Status: Post processing for IpsoOS ...  
>> Execute cmd: LAUNCH=</bin/bash /tftpboot/install-services >  
>> Execute cmd: options log=0 canfail=0 cont=0 argc=2>  
>> Execute cmd: pwd=</tftpboot>
```

```
*** ERROR: ***
```

Could not enable service tftp: not in xinetd.conf

Failed to get unit file state for dhcpd.service: No such file or directory

Failed to execute operation: No such file or directory

(systemctl enable dhcpd.service) [FAILED]

sed: can't read /usr/lib/systemd/system/dhcpd.service: No such file or directory

/tftpboot/service-lib.sh: line 1272: restorecon: command not found

\*\*\* ERROR: \*\*\*

restorecon -v /etc/systemd/system/dhcpd.service

/tftpboot/service-lib.sh: line 1272: restorecon: command not found

\*\*\* ERROR: \*\*\*

restorecon -v /etc/dhcpd.conf

Failed to start dhcpd.service: Unit dhcpd.service failed to load: Invalid argument. See system logs and 'systemctl status dhcpd.service' for details.

(systemctl start dhcpd.service) [FAILED]

(systemctl status dhcpd.service) [FAILED]

Output:

Note: Using init directory /etc/init.d

Note: Using Systemd for service control

\*\*\* WARNING: \*\*\*

network interface eth1 not found

/tftpboot/install-services: install services: tftp dhcpd

systemctl enable dhcpd.service ... command: echo > /tftpboot/dhcpd.service

command: cp -af /tftpboot/dhcpd.service /etc/systemd/system/dhcpd.service

command: restorecon -v /etc/systemd/system/dhcpd.service

command: cp -af /tftpboot/dhcpd.conf /etc/dhcpd.conf

command: restorecon -v /etc/dhcpd.conf

systemctl -q daemon-reload ... [OK]

systemctl start dhcpd.service ... systemctl status dhcpd.service ... \*

dhcpd.service

Loaded: error (Reason: Invalid argument)

Active: inactive (dead)

May 18 11:51:47 cie-55-31 systemd[1]: dhcpd.service lacks both ExecStart= and ExecStop= setting. Refusing.

May 18 11:51:47 cie-55-31 systemd[1]: dhcpd.service lacks both ExecStart= and ExecStop= setting. Refusing.

/tftpboot/install-services: terminated with error status 1

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```
during command /tftpboot/install-services
Errors:
/tftpboot/service-lib.sh: line 241: rpm: command not found
*** Fatal error: ***

/tftpboot/install-services: is not a number

Output:
/tftpboot/install-services: terminated with status 99
```

i

```
Error: Execute cmd: exit=99
Error: Execute cmd: error=</usr/diskless/crco_data/service-lib.sh: line 241:
rpm: command not found
*** Fatal error: ***

/usr/diskless/crco_data/crco-install: is not a number
>
Error: Execute cmd: exit=99
Error: Execute cmd: error=</tftpboot/service-lib.sh: line 241: rpm: command
not found
*** Fatal error: ***

/tftpboot/install-services: is not a number
```

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```
during command installation script (post)
Errors:
./Bruker/service-lib.sh: line 241: rpm: command not found
is not a number
/home/sermnadmin/Downloads/Topspin-3.2.7/Topspin3.2PL7/xwinstall.d/modscript
s/flexlm.post: terminated with status 99
```

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```
*** Error ***
*
Could not show firewall rules for Diskless:
(command: /bin/bash /usr/diskless/prog/bin/setfirewall.sh show)

/usr/diskless/prog/bin/service-lib.sh: line 241: rpm: command not found
Fatal error: is not a number
/usr/diskless/prog/bin/setfirewall.sh: terminated with status 3
```

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(arquitectura i386)

```
sermnadmin@cie-55-31:~$ topspin
TopSpin 3.5 pl 6 - Copyright (C) 2016 Bruker BioSpin GmbH
Installation directory: "/opt/topspin3.5pl6"
Waiting for FLEXlm license
```

```
Cannot checkout TopSpin FLEXlm license
Invalid license file syntax
Feature:      TOPSPIN3
License path: /usr/local/flexlm/Bruker/licenses/license.dat
FLEXlm error: -2,413
For further information, refer to the FLEXlm End User Manual,
available at "www.macrovision.com".
```

The FLEXlm host ID of this machine is 0010182755de

If your workstation controls the spectrometer,  
you may run the program without license.  
This requires the acquisition server to be ready.

checking connection to spectrometer (may last several minutes)

```
CPR : 2017-05-19 17:30:51.941 +0200
TopSpin is currently running under
the spectrometer service license.
Please apply to receive a correct
license for your workstation
using the online license form:
https://www.bruker.com/nmr_license_requests.html
```

```
error message cprserver: GetSystemPrinterList: dlopen(libcups.so) failed
Exception in thread "main" java.lang.UnsatisfiedLinkError:
/opt/topspin3.5pl6/jre/lib/i386/libawt_xawt.so: libXtst.so.6: cannot open
shared object file: No such file or directory
```

```
at java.lang.ClassLoader$NativeLibrary.load(Native Method)
at java.lang.ClassLoader.loadLibrary0(ClassLoader.java:1941)
at java.lang.ClassLoader.loadLibrary(ClassLoader.java:1824)
at java.lang.Runtime.load0(Runtime.java:809)
at java.lang.System.load(System.java:1086)
at java.lang.ClassLoader$NativeLibrary.load(Native Method)
at java.lang.ClassLoader.loadLibrary0(ClassLoader.java:1941)
at java.lang.ClassLoader.loadLibrary(ClassLoader.java:1845)
at java.lang.Runtime.loadLibrary0(Runtime.java:870)
at java.lang.System.loadLibrary(System.java:1122)
at java.awt.Toolkit$3.run(Toolkit.java:1636)
at java.awt.Toolkit$3.run(Toolkit.java:1634)
at java.security.AccessController.doPrivileged(Native Method)
at java.awt.Toolkit.loadLibraries(Toolkit.java:1633)
at java.awt.Toolkit.<clinit>(Toolkit.java:1668)
at java.awt.Color.<clinit>(Color.java:275)
at de.bruker.nmr.mfw.base.BSplashScreen.<clinit>(BSplashScreen.java:44)
```

```
at de.bruker.nmr.mfw.base.MfwBuilder.programMain(MfwBuilder.java:420)
at bruker.bio.start.topspin.Start.main(Start.java:47)
```

premature Java Virtual Machine termination  
Program is exiting ...

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Ara el TopSpin ja s'executa,

```

sermnadmin@cie-55-31:~$ topspin
TopSpin 3.5 pl 6 - Copyright (C) 2016 Bruker BioSpin GmbH
Installation directory: "/opt/topspin3.5pl6"
Waiting for FLEXlm license

Cannot checkout TopSpin FLEXlm license
Invalid license file syntax
Feature:      TOPSPIN3
License path: /usr/local/flexlm/Bruker/licenses/license.dat
FLEXlm error: -2,413
For further information, refer to the FLEXlm End User Manual,
available at "www.macrovision.com".

The FLEXlm host ID of this machine is 0010182755de

If your workstation controls the spectrometer,
you may run the program without license.
This requires the acquisition server to be ready.

checking connection to spectrometer (may last several minutes)
CPR : 2017-05-19 17:33:57.165 +0200
TopSpin is currently running under
the spectrometer service license.

```

Please apply to receive a correct  
license for your workstation  
using the online license form:  
[https://www.bruker.com/nmr\\_license\\_requests.html](https://www.bruker.com/nmr_license_requests.html)

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IPSO: connected to spectrometer subnet

- TCP/IP address = 149.236.99.90
- IPSO type: AQS
- Tctrl : 1
- Fctrls: 2
- Gctrl1: with digital preemphasis
- Rctrl : 1
- FREDs : none
- DPP : 1

DRU1: AQS DRU Z100977/00582 ECL 04.00

- TCP/IP address = 149.236.99.89
- Firmware Version = 151014
- DRU1 controls AQS-Rack and HPPR/2

AQS: connected to 149.236.99.89:/dev/tty10

\_Slot\_ SBSB \_\_\_\_\_ Board \_\_\_\_\_

| Number | Addr | Type | HW-VS | FW-VS | ID | ECL | Name   | Description                                      |
|--------|------|------|-------|-------|----|-----|--------|--------------------------------------------------|
| 2      | 0x10 | 0x72 | 0x3   | AW    | R  | 3.0 | REC-1  | AQS RXAD400 Z102116/705 ECL                      |
| 3      | 0x34 | 0x2  | 0x1   |       | X  | 5.0 | REF-1  | REF-400 Reference Board for AQS Receiver         |
| 4      | 0x24 | 0x11 | 0x8   | AE    | S  | 2.0 | SGU-1  | AQS SGU/2 400 Z103080/00351 ECL 02.00 FPGA-Vs=AF |
| 5      | 0x25 | 0x11 | 0x8   | AE    | S  | 2.0 | SGU-2  | AQS SGU/2 400 Z103080/00340 ECL 02.00 FPGA-Vs=AF |
| 9      | 0x36 | 0x5  | 0     |       | A  | 6.0 | BLA-1  | BLA2BB 150/60 20-400                             |
| 21     | 0x21 | 0xcf | 0     |       | P  | 1.0 | PSD-1  | PSD/3 Power Supply Distribution Board            |
| --     | 0x20 | 0xd1 | 0     |       | B  | 2.0 | MASTER | AQS/3 Chassis                                    |
| 1      | 0x20 | 0x7  | 0     |       | B  |     | MASTER | AQS Rack Master                                  |

Logical Receiver/SGU configuration:

Router: 1 AQS-Minirouter

Amplifiers (AQS configuration):

| R0 | Amplifier | Module | Nucleus | Power/W | Switchbox | Name         | PN/SN/FW      |
|----|-----------|--------|---------|---------|-----------|--------------|---------------|
| 1  | 1         | 1      | B       | 150     | N         | BLA2BB       | 150/60 20-400 |
| 2  | 1         | 2      | B       | 60      | N         | BLA2BB       | 150/60 20-400 |
| 3  | 2         | 1      | 2H      | 20      | N         | 2HTX in BSMS | -/-/-         |

Gradient amplifiers at the spectrometer subnet:

BGA1: BGA\_W1213891\_0040

- TCP/IP address = 149.236.99.93
- Firmware version = 20080206
- Web version = 1.1
- Current limits = 0.0/X, 0.0/Y, -10.0/Z (in A)

BSMS: BSMS/2 connected to ethernet

- TCP/IP address = 149.236.99.20
- ELCB firmware version = 20150423
- SCB channels = 20
- Shim System = BOSS1-SB
- 2HTX Amplifier: available
- Lock: installed

MAS2 control unit: connected to 149.236.99.90:/dev/tty05

- Boot Software : 20031216
- Application Software : 20120329
- Insert Time : 10 s
- Eject Time : 10 s
- Spin Lock Tolerance : 5 Hz

Eurotherm VTU: device connected to 149.236.99.90:/dev/tty04

Preamplifiers :

HPPR2: - HPPR/2 preamplifier connected to 149.236.99.89:/dev/tty10  
Type : HPPR/2  
Controller: Cover/1 with Application firmware = AQ  
15 LEDs for tuning, 15 LEDs for matching  
Module 1 : 1H/2H 1H (virtual 50 Ohm reference: 274.9%/85.9deg)  
PN=Z003959, SN=00418 from 20070319  
Module 2 : 1H/2H 2H (virtual 50 Ohm reference: 116.6%/-71.1deg)  
PN=Z003959, SN=00418 from 20070319  
Module 3 : XBB19F 2HS  
PN=Z003504, SN=00415 from 20070425

Gradient Controller cable connections

-----  
GCtrl1: connected via DPP to GREAT 1-10 Gradient Amplifier Unit  
W1213891/0040

RF cable connections (assumed)

-----  
SGU1 NORM output -> input 2 of transmitter 1 (BLA2BB 150/60 20-400  
W1345049/1817)  
SGU1 AUX output -> input 1 of transmitter 2 (2HTX in BSMS)  
SGU2 NORM output -> input 3 of transmitter 1 (BLA2BB 150/60 20-400  
W1345049/1817)  
SGU2 AUX output -> TUNE signal input of HPPR

Blanking cable connections (assumed)

-----  
transmitter 1 = BLA2BB 150/60 20-400 W1345049/1817:  
- amplifier B-150W uses blanking 1  
- amplifier B-60W uses blanking 2  
  
transmitter 2 = 2HTX in BSMS:  
- amplifier 2H-20W uses blanking 3

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```
atma_ProbeCorba_impl::start exception: Error:can't stat the wobble process  
[NucBB-NucBB-]
```

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```
Atma Info  
atma_ProbeCorba_impl::start exception: Error:can't start the wobble process  
[NucBB-NucBB-]  
(Original message = Atma Info  
atma_ProbeCorba_impl::start exception: Error:can't start the wobble process  
[NucBB-NucBB-])  
  
=====  
atma_ProbeCorba_impl::start exception: Error:can't start the wobble process  
[NucBB-NucBB-]  
  
24 May 2017 11:30:08.978 +0200  
Topspin Version  = 3.5 pl 6 (of August 4 2016),build 1784  
JVM Version      = 1.8.0_102 (32 bit) Oracle Corporation  
JVM Total memory = 342 MB  
JVM Free  memory = 189 MB  
  
java.lang.Exception: Stack trace  
    at  
bruker.bio.topspin.bfw.gui.warn.AbstractMessage.getDetailsString(AbstractMes  
sage.java:827)  
    at  
bruker.bio.topspin.bfw.gui.warn.AbstractMessage.setMsgSource(AbstractMessage  
.java:790)  
    at  
bruker.bio.topspin.bfw.gui.warn.AbstractMessage.<init>(AbstractMessage.java:  
774)  
    at de.bruker.nmr.mfw.base.BWarning.<init>(BWarning.java:102)  
    at de.bruker.nmr.mfw.base.BError.<init>(BError.java:93)  
    at  
de.bruker.nmr.sc.atma.AtmaPanelStart.showAtmaException(AtmaPanelStart.java:1  
972)  
    at de.bruker.nmr.sc.atma.AtmaPanelStart.connect(AtmaPanelStart.java:280)
```

```
at de.bruker.nmr.sc.atma.AtmaPanelStart.start(AtmaPanelStart.java:138)
at de.bruker.nmr.sc.atma.AtmaPanelStart.atmm(AtmaPanelStart.java:110)
at sun.reflect.NativeMethodAccessorImpl.invoke0(Native Method)
at
sun.reflect.NativeMethodAccessorImpl.invoke(NativeMethodAccessorImpl.java:62)
at
sun.reflect.DelegatingMethodAccessorImpl.invoke(DelegatingMethodAccessorImpl.java:43)
at java.lang.reflect.Method.invoke(Method.java:498)
at
de.bruker.nmr.mfw.base.CmdThreadImpl.doJmethod(CmdThreadImpl.java:600)
at de.bruker.nmr.mfw.base.CmdThreadImpl.exec(CmdThreadImpl.java:509)
at de.bruker.nmr.mfw.base.CmdThreadImpl.run(CmdThreadImpl.java:489)
```

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- High Resolution Systems
- Solid State Systems
- Micro-Imaging and Diffusion Systems

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```
nmr_save.c:30:19: fatal error:stdio.h: No such file or directory
#include <stdio.h>
                ^
compilation terminated
```

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|                                   |                                                                                     |
|-----------------------------------|-------------------------------------------------------------------------------------|
| <b>Customer Info</b>              |                                                                                     |
| <b>Postal Code</b>                | 08193                                                                               |
| <b>City *</b>                     | Cerdanyola del Vallès                                                               |
| <b>Country *</b>                  | Catalunya (Spain)                                                                   |
| <b>Phone Contact Customer *</b>   | +34 93 581 2291                                                                     |
| <b>Mobile</b>                     | —                                                                                   |
| <b>Fax</b>                        | —                                                                                   |
| <b>E-Mail *</b>                   | <a href="mailto:s.rmn@uab.cat">s.rmn@uab.cat</a>                                    |
| <b>Bruker Info</b>                |                                                                                     |
| <b>Engineer *</b>                 | Òscar Benítez                                                                       |
| <b>Office *</b>                   | Bruker Española, S.A.                                                               |
| <b>Central Hotline Phone *</b>    | +34 686 944 920                                                                     |
| <b>Central Hotline E-Mail *</b>   | <a href="mailto:oscar.benitez@bruker.com">oscar.benitez@bruker.com</a>              |
| <b>System Info</b>                |                                                                                     |
| <b>Order No. *</b>                |    |
| <b>Contract Service No</b>        | —                                                                                   |
| <b>System Type</b>                | Avance III 400                                                                      |
| <b>Console Part and Serial No</b> |    |
| <b>Coil</b>                       |    |
| <b>Dewar</b>                      |   |
| <b>Shim System Offset *</b>       |  |
| <b>Shim System Angle</b>          | —                                                                                   |
| <b>CryoProbe Order No.</b>        | —                                                                                   |
| <b>Location</b>                   | Servei de Ressonància Magnètica Nuclear                                             |
| <b>Register No.</b>               |  |

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-  <TOPSPIN\_HOME>/exp/stan/nmr/par/users/
-  <TOPSPIN\_HOME>/exp/stan/nmr/par/teo\* (i pau\*, mir\*. Tots es copien a <TOPSPIN\_HOME>/exp/stan/nmr/par/users/)
-  <TOPSPIN\_HOME>/exp/stan/nmr/lists/pp/users/
-  <TOPSPIN\_HOME>/exp/stan/nmr/lists/gp/users/ (carpeta buida)
-  <TOPSPIN\_HOME>/exp/stan/nmr/lists/vc/ (carpeta buida)
-  <TOPSPIN\_HOME>/exp/stan/nmr/lists/vd/
-  <TOPSPIN\_HOME>/exp/stan/nmr/lists/vt/ (carpeta buida)
-  <TOPSPIN\_HOME>/exp/stan/nmr/lists/wave/users/

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```
# 20160422 - MCE - Hide nmrsu user from the login screen
IncludeAll = false
Include = sermnadmin,sermnuab
```

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```
# 20160422 - MCE - Prevent nmrsu user from log-in via ssh
DenyUsers nmrsu
```

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```
root@cie-55-31:/home/sermnadmin/Documents/old-old-hdd-recovery/sdb4-old-old-hdd-opt/topspin/data# rsync -Hvax --progress --stats ./*/
/opt/topspin3.5pl6/data/
[...]
Number of files: 7,573 (reg: 6,405, dir: 1,168)
Number of created files: 7,573 (reg: 6,405, dir: 1,168)
Number of deleted files: 0
Number of regular files transferred: 6,405
Total file size: 459,101,587 bytes
Total transferred file size: 459,101,587 bytes
Literal data: 459,101,587 bytes
Matched data: 0 bytes
File list size: 65,536
File list generation time: 0.001 seconds
File list transfer time: 0.000 seconds
Total bytes sent: 459,608,752
Total bytes received: 128,851

sent 459,608,752 bytes  received 128,851 bytes  54,086,776.82 bytes/sec
total size is 459,101,587  speedup is 1.00
```

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**WORK-IN-PROGRESS** Started on Monday 22 May, at 16:41

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```
root@cie-55-31:/home/sermnadmin/Documents/old-old-hdd-recovery/sdb4-old-old-hdd-opt/sermn-uab# rsync -Hvax --progress --stats ./data /home/sermnuab/
```

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```
[root@cie-55-32 etc]# cat /etc/exports
# export data folder to SeRMN data server
/home/sermnuab/data
sermn.uab.es(ro,sync,no_subtree_check,root_squash,all_squash,insecure)
#
# export lists folder to SeRMN data server
/opt/topspin3.2/exp/stan/nmr/lists
sermn.uab.es(ro,sync,no_subtree_check,root_squash,all_squash,insecure)
[...]
data/test/nmr/tune-31P/1/pdata/1/title
13 100%    0.02kB/s    0:00:00 (xfr#1048377, to-chk=0/1245348)

Number of files: 1,245,348 (reg: 1,048,377, dir: 196,971)
Number of created files: 1,245,348 (reg: 1,048,377, dir: 196,971)
Number of deleted files: 0
Number of regular files transferred: 1,048,377
Total file size: 85,823,932,533 bytes
Total transferred file size: 85,823,932,533 bytes
Literal data: 85,823,932,533 bytes
Matched data: 0 bytes
File list size: 4,915,041
File list generation time: 0.001 seconds
File list transfer time: 0.000 seconds
Total bytes sent: 85,909,762,834
Total bytes received: 21,195,209

sent 85,909,762,834 bytes   received 21,195,209 bytes   23,906,233.98
bytes/sec
total size is 85,823,932,533   speedup is 1.00
```

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|        | Probe: 1H-BB/19F BBO+ATMA     |                |                    | Console: SGU & Amp         |
|--------|-------------------------------|----------------|--------------------|----------------------------|
| Nuclei | ATMA / ATMM<br>(Tune & Match) | Power<br>Limit | PROSOL<br>(Pulses) | CORTAB<br>(RF Calibration) |
| 1H     | ✓                             | ✓              | ✓                  | ✓                          |
| 13C    | ✓                             | ✓              | ✓                  | ✓                          |
| 31P    | ✓                             | ✓              | ✓                  | ✓                          |

From:

<https://sermn.uab.cat/wiki/> - SeRMN Wiki

Permanent link:

[https://sermn.uab.cat/wiki/doku.php?id=informatica:hp\\_xw4400\\_400sb\\_topspin35&rev=1496141701](https://sermn.uab.cat/wiki/doku.php?id=informatica:hp_xw4400_400sb_topspin35&rev=1496141701)

Last update: 2017/05/30 12:55