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```
sermnadmin@cie-55-32:~/Downloads/Topspin-3.2.7/Topspin3.2PL7$ ./install.cmd
NOTE: using DISPLAY=:0
Please enter root password if prompted for it.
Password:
NOTE: using DISPLAY=:0
NOTE: using
TCL_LIBRARY=/home/sermnadmin/Downloads/Topspin-3.2.7/Topspin3.2PL7/linux/tcl-8.5.11/lib/tcl8.5
NOTE: using
TK_LIBRARY=/home/sermnadmin/Downloads/Topspin-3.2.7/Topspin3.2PL7/linux/tk-8.5.11/lib/tk8.5
NOTE: running
/home/sermnadmin/Downloads/Topspin-3.2.7/Topspin3.2PL7/linux/tk-8.5.11/bin/wish8.5 -f
/home/sermnadmin/Downloads/Topspin-3.2.7/Topspin3.2PL7/xwinstall.d/swim/lib/xwinstall.tcl --
/home/sermnadmin/Downloads/Topspin-3.2.7/Topspin3.2PL7/linux/tk-8.5.11/bin/wish8.5: error while loading shared libraries: libX11.so.6: cannot open shared object file: No such file or directory
Installation program failed with status 127
(Command:
/home/sermnadmin/Downloads/Topspin-3.2.7/Topspin3.2PL7/linux/tk-8.5.11/bin/wish8.5 -f
/home/sermnadmin/Downloads/Topspin-3.2.7/Topspin3.2PL7/xwinstall.d/swim/lib/xwinstall.tcl --)
```

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```
during command /tftpboot/install-services
Errors:
/tftpboot/service-lib.sh: line 241: rpm: command not found
*** Fatal error: ***

/tftpboot/install-services: is not a number

Output:
/tftpboot/install-services: terminated with status 99
```

i

```
Error: Execute cmd: exit=99
Error: Execute cmd: error=</usr/diskless/crco_data/service-lib.sh: line 241:
rpm: command not found
*** Fatal error: ***

/usr/diskless/crco_data/crco-install: is not a number
>
Error: Execute cmd: exit=99
Error: Execute cmd: error=</tftpboot/service-lib.sh: line 241: rpm: command
not found
*** Fatal error: ***
```

```
/tftpboot/install-services: is not a number
```

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```
during command installation script (post)
Errors:
./Bruker/service-lib.sh: line 241: rpm: command not found
is not a number
/home/sermnadmin/Downloads/Topspin-3.2.7/Topspin3.2PL7/xwinstall.d/modscript
s/flexlm.post: terminated with status 99
```

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```
*** Error ***
*
Could not show firewall rules for Diskless:
(command: /bin/bash /usr/diskless/prog/bin/setfirewall.sh show)

/usr/diskless/prog/bin/service-lib.sh: line 241: rpm: command not found
Fatal error: is not a number
```

```
/usr/diskless/prog/bin/setfirewall.sh: terminated with status 3
```

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```
mount swapfile ASP_ST2:/usr/diskless/clients/spect/swapfile failed, rpc
status 2 RPC: Can't decode result
```

```
PANIC: nfs_swapvp failed mounting swap or dump device
```

```
syncing disks... done
```

```
Fast ethernet (FEN): 100 MBit/s, half duplex
```

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```
RISC/os Release 4_52 mips Version UMIPS
```

```
Total real memory = 16777216
Available memory   = 14442496
Root on nfs file :, Swap on nfs file :
Fast ethernet (FEN) : 100 MBit/s, full duplex
hostname: spect
domainname: (none)
Root fstype nfs
Available memory   = 12795904

Fast ethernet (FEN) : 100 MBit/s, full duplex
```

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```
Running Power-On Diagnostics...
Low Memory Test... PASSED
ID PROM Test... PASSED
Write Buffer Test... PASSED
Memory Tests... PASSED
SCC Test... PASSED
Timer/Time-of-Day Clock Test... *** SKIPPED ***
Battery Check Test...state 00000002
PASSED
NVRAM Test... PASSED
Lance Register Test ... *** SKIPPED ***
Lance Master Test ... *** SKIPPED ***

Fast ethernet (FEN): 100 MBit/s, half duplex

Power-On Diagnostics..... PASSED
  Main Memory Test..... PASSED  ID PROM Test.....
PASSED
  Write Buffer Test..... PASSED  SCC Tests.....
PASSED
  NVRAM Tests..... PASSED  Ethernet Tests.....
PASSED

BRUKER CCU R4600 Monitor
Version 5.40 MIPS OPT Thu May 17 12:18:17 MSZ 2001 mar
Memory size: 16777216 (0x1000000) bytes, 16 MB
Icache size: 16384 (0x4000) bytes
```

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```
Apr 21 15:22:17 INIT: Command is respawning too rapidly. Check for possible errors.  
id: aq "/etc/startd"
```

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```
spect > more /etc/inittab  
...  
aq:234:respawn:/etc/startd  
ec:234:off:/etc/ecud  
...
```

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```
sermadmin@cie-55-32:/opt/topspin3.2$ ./topspin  
TopSpin 3.2 pl 7 - Copyright (C) 2015 Bruker BioSpin GmbH  
Installation directory: "/opt/topspin3.2"  
Waiting for FLEXlm license
```

```
2016-04-21T13:59:59.900+02:00
Cannot checkout TopSpin FLEXlm license
FEATURE TOPSPIN3 Invalid license file syntax
Feature:          TOPSPIN3
License path:    /usr/local/flexlm/Bruker/licenses/license.dat
FLEXlm error:   -2,413
For further information, refer to the FLEXlm End User Manual,
available at "www.macrovision.com".
```

The FLEXlm host ID of this machine is a0369f8f0ab3

If your workstation controls the spectrometer,
you may run the program without license.
This requires the acquisition server to be ready.

```
checking connection to spectrometer (may last a minute)
no spectrometer service license available
```

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```
# SerMN - Servers and network devices
#
[...]
158.109.56.84    rmn6.uab.cat      rmn6              # license server
```

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```
sermnadmin@cie-55-32:/opt/topspin3.2$ topspin
TopSpin 3.2 pl 7 - Copyright (C) 2015 Bruker BioSpin GmbH
Installation directory: "/opt/topspin3.2"
Waiting for FLEXlm license

2016-04-21T16:53:36.534+02:00
Cannot checkout TopSpin FLEXlm license
FEATURE TOPSPIN3 Invalid license file syntax
Feature:          TOPSPIN3
License path:    /usr/local/flexlm/Bruker/licenses/license.dat
FLEXlm error:   -2,413
For further information, refer to the FLEXlm End User Manual,
available at "www.macrovision.com".
```

The FLEXlm host ID of this machine is a0369f8f0ab3

If your workstation controls the spectrometer,
you may run the program without license.
This requires the acquisition server to be ready.

```
checking connection to spectrometer (may last a minute)
CPR : 2016-04-21 16:53:36.758 +0200
  TopSpin is currently running under
  the spectrometer service license.
  Please apply to receive a correct
  license for your workstation
  using the online license form:
  https://www.bruker.com/nmr_license_requests.html
```

```
error message cprserver: GetSystemPrinterList: dlopen(libcups.so) failed
Exception in thread "main" java.lang.UnsatisfiedLinkError:
/opt/topspin3.2/jre/lib/i386/libawt_xawt.so: libXext.so.6: cannot open
shared object file: No such file or directory
  at java.lang.ClassLoader$NativeLibrary.load(Native Method)
  at java.lang.ClassLoader.loadLibrary0(ClassLoader.java:1937)
  at java.lang.ClassLoader.loadLibrary(ClassLoader.java:1822)
  at java.lang.Runtime.load0(Runtime.java:809)
  at java.lang.System.load(System.java:1083)
  at java.lang.ClassLoader$NativeLibrary.load(Native Method)
  at java.lang.ClassLoader.loadLibrary0(ClassLoader.java:1937)
  at java.lang.ClassLoader.loadLibrary(ClassLoader.java:1843)
  at java.lang.Runtime.loadLibrary0(Runtime.java:870)
  at java.lang.System.loadLibrary(System.java:1119)
  at java.awt.Toolkit$3.run(Toolkit.java:1636)
  at java.awt.Toolkit$3.run(Toolkit.java:1634)
  at java.security.AccessController.doPrivileged(Native Method)
  at java.awt.Toolkit.loadLibraries(Toolkit.java:1633)
  at java.awt.Toolkit.<clinit>(Toolkit.java:1668)
  at java.awt.Color.<clinit>(Color.java:275)
  at de.bruker.nmr.mfw.root.FramePars.<init>(FramePars.java:82)
  at de.bruker.nmr.mfw.base.Start.<clinit>(Start.java:55)
```

```
premature Java Virtual Machine termination
Program is exiting ...
```

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Ara el TopSpin ja s'executa,

```
sermnadmin@cie-55-32:/opt/topspin3.2$ topspin
TopSpin 3.2 pl 7 - Copyright (C) 2015 Bruker BioSpin GmbH
Installation directory: "/opt/topspin3.2"
Waiting for FLEXlm license

2016-04-22T10:27:59.081+02:00
Cannot checkout TopSpin FLEXlm license
FEATURE TOPSPIN3 Invalid license file syntax
Feature:          TOPSPIN3
License path:    /usr/local/flexlm/Bruker/licenses/license.dat
FLEXlm error:   -2,413
For further information, refer to the FLEXlm End User Manual,
available at "www.macrovision.com".

The FLEXlm host ID of this machine is a0369f8f0ab3

If your workstation controls the spectrometer,
you may run the program without license.
This requires the acquisition server to be ready.

checking connection to spectrometer (may last a minute)
```

```
CPR : 2016-04-22 10:27:59.296 +0200
TopSpin is currently running under
the spectrometer service license.
Please apply to receive a correct
license for your workstation
using the online license form:
https://www.bruker.com/nmr_license_requests.html
```

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Amplifier	Preamplifier	Probe: 4 mm MAS 1H/BB H8113/0028 [33]
[X 100 W]	[HPHP XBB31P]	[BB] 2 Inner Coil
[H 50 W]	[HPHP 19F/1H]	[1H] 1 Inner Coil

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- High Resolution Systems
- Solid State Systems
- Micro-Imaging and Diffusion Systems

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```
Wrong version of prosol files! Call edprosol to update it!  
  
You may start a read only session of the editor in order to save your own  
modifications.  
  
Or you may start an automatic update of your prosol directory branch,  
wherein all prosol parameter files are replaced. Your old 90 deg. pulses are  
saved and the new pulse files are recalculated based on them.
```

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- Read only editor for individual saving
- Automatic Update, saving only 90 deg. reference pulses.

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```
# 20160422 - MCE - Hide nmrsu user from the login screen
IncludeAll = false
Include = sermnadmin,sermnub
```

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```
# 20160422 - MCE - Prevent nmrsu user from log-in via ssh
DenyUsers nmrsu
```

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```
[root@cie-55-32 etc]# cat /etc/exports
# export data folder to SeRMN data server
/home/sermnuab/data
sermn.uab.es(ro, sync, no_subtree_check, root_squash, all_squash, insecure)
#
# export lists folder to SeRMN data server
/opt/topspin3.2/exp/stan/nmr/lists
sermn.uab.es(ro, sync, no_subtree_check, root_squash, all_squash, insecure)

[...]
```

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```
nmr_save.c:30:19: fatal error:stdio.h: No such file or directory
#include <stdio.h>
                ^
compilation terminated
```

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solució al 600.

L'error és:

```
ILLEGAL TERMINATION
of command /opt/topspin3.2/prog/mod/xwish3
MODULE DIED IN ITS SLEEP
due to signal 11 (Segmentation fault)
(Original message = ILLEGAL TERMINATION
of command /opt/topspin3.2/prog/mod/xwish3
MODULE DIED IN ITS SLEEP
due to signal 11 (Segmentation fault))

=====
14 August 2019 13:00:44.137 +0200
Topspin Version = 3.2 (of July 20 2015),build 1791
JVM Version      = 1.8.0_31 Oracle Corporation
JVM Total memory = 174 MB
JVM Free  memory = 121 MB

java.lang.Exception: Stack trace
    at
de.bruker.nmr.mfw.base.AbstractMessage.setMsgSource(AbstractMessage.java:813
)
    at
de.bruker.nmr.mfw.base.AbstractMessage.<init>(AbstractMessage.java:772)
    at de.bruker.nmr.mfw.base.BWarning.<init>(BWarning.java:57)
    at
de.bruker.nmr.prsc.cpr.CprListener$CPRWarning.<init>(CprListener.java:337)
    at de.bruker.nmr.prsc.cpr.CprListener.loadAlert(CprListener.java:878)
    at de.bruker.cpr.CPRClientPOA._invoke_loadAlert(CPRClientPOA.java:121)
    at de.bruker.cpr.CPRClientPOA._invoke(CPRClientPOA.java:43)
    at org.openorb.adapter.poa.POA.dispatch(Unknown Source)
    at org.openorb.net.AbstractServerRequest.dispatch(Unknown Source)
    at org.openorb.net.ServerManagerImpl.serve_request(Unknown Source)
    at org.openorb.net.ServerManagerImpl.thread_pool_main(Unknown Source)
    at org.openorb.net.ServerManagerImpl.access$200(Unknown Source)
    at org.openorb.net.ServerManagerImpl$PoolThread.run(Unknown Source)
```

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