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- Select the type of acquisition: High Resolution Systems
- Available spectrometer or datastation configurations
  - Select the configuration you want to use: spect
- Select the items you want to install
  - Install Pulse Programs
  - Install Bruker AU Programs
  - [unchecked] Recompile All User AU Programs
  - Install Library CPD Programs
  - Install Library Gradient Files
  - Install Library Shape Files
  - Convert Standard Parameter Sets
  - Install Standard Scaling Region Files
  - Install Bruker Python Programs
- Select the basic frequency of your spectrometer
  - Basic Frequency: 600.13
- Select the pre-scan delay DE
  - Default pre-scan-delay (us): 6.5
- Select the plotter paper format
  - Paper format: A4 / Letter

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- SAVE INSTALLATION FILES
  - Location of backup file: /opt/topspin3.6.3/nmr\_backup \* Installation to be saved (TopSpin home): /opt/topspin3.6.3
  - Spectrometer configuration (e.g. spect): spect
  - Display default informaiton: checked
  - Display additional informaiton: unchecked
- Execute "Save installation specific files" periodically: [Automatic Backup]

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```
retrieving path to diskless (may take some time)...
found /usr/diskless
prepare backup: Copy /usr/diskless/prog/logfiles to
/opt/topspin3.6.3/nmr_backup/tmp_backup_dir/logfiles
extract and backup /opt/topspin3.6.3/conf/instr/servtool/bsmstool/boss/
```

```
=>
extract and backup /opt/topspin3.6.3/conf/instr/servtool/bsmstool/boss/wb/
now extracting from dir
/opt/topspin3.6.3/conf/instr/servtool/bsmstool/boss/wb/...

<=
=>
extract and backup
/opt/topspin3.6.3/conf/instr/servtool/bsmstool/boss/boss3_sb/
now extracting from dir
/opt/topspin3.6.3/conf/instr/servtool/bsmstool/boss/boss3_sb/...

<=
now extracting from dir
/opt/topspin3.6.3/conf/instr/servtool/bsmstool/boss/...

extract and backup /opt/topspin3.6.3/exp/stan/nmr/py/
=>
extract and backup /opt/topspin3.6.3/exp/stan/nmr/py/biotop/
=>
extract and backup /opt/topspin3.6.3/exp/stan/nmr/py/biotop/experiments/
[...]...
now extracting from dir /opt/topspin3.6.3/plot/layouts/layouts.stack/...
now extracting from dir /opt/topspin3.6.3/exp/stan/nmr/lists/scl/...

Now writing zip file
/opt/topspin3.6.3/nmr_backup/extractedDirectories.zip1...

Every file of the following directories will be saved:

/opt/topspin3.6.3/QTP/
/opt/topspin3.6.3/conf/global/
/opt/topspin3.6.3/conf/instr/...
[...]
/opt/topspin3.6.3/exp/stan/nmr/lists/ ...
[...]
/opt/topspin3.6.3/db/quantref/
/opt/topspin3.6.3/classes/prop/ ...
[...]
/opt/topspin3.6.3/db/user/
/opt/topspin3.6.3/data/final/nmr/protocolfiles/

Additionally the following files will be saved (if existing):

/opt/topspin3.6.3/conf/instr/spect/uxnmr.par
/opt/topspin3.6.3/conf/instr/spect/uxnmr.info
[...]
/opt/topspin3.6.3/conf/instr/spect/hardware_list
/opt/topspin3.6.3/classes/prop/ ...
[...]
/opt/topspin3.6.3/conf/instr/ ...
```

```
[...]  
/opt/topspin3.6.3/conf/topspin-users.prop  
/opt/topspin3.6.3/exp/stan/nmr/lists/probeheads  
/opt/topspin3.6.3/prog/server/export.conf  
  
-----  
  
start to generate back-up copy...  
  
Now writing zip file  
/opt/topspin3.6.3/nmr_backup/completeDirectories.zip1...  
  
Now writing zip file  
/opt/topspin3.6.3/nmr_backup/completeDirectories.zip2...  
  
Now writing zip file  
/opt/topspin3.6.3/nmr_backup/nmr_backup_20210812-1054.zip...  
  
#####  
  
Backing up of NMR data was successful!  
The backup file  
  
/opt/topspin3.6.3/nmr_backup/nmr_backup_20210812-1054.zip  
  
has been generated!  
  
#####
```

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- SAVE USER FILES
  - Location of backup file: /opt/topspin3.6.3/nmr\_backup
  - Overwrite existing backup file unchecked
  - Display default information: checked
  - Display additional information: unchecked
- Execute Save user files periodically:

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```
retrieving path to diskless (may take some time)...
```

found /usr/diskless

Every non-Bruker file in /home/sermnadmin/.topspin-cie-56-18.uab.cat/ will be saved:

Every non-Bruker file in /opt/topspin3.6.3/exp/stan/nmr/au/src/ will be saved:

now extracting from dir /opt/topspin3.6.3/exp/stan/nmr/au/src/...

Now writing zip file

/opt/topspin3.6.3/nmr\_backup/extractedDirectories.zip1...

Every non-Bruker file in /opt/topspin3.6.3/exp/stan/nmr/lists/base\_info/ will be saved:

now extracting from dir /opt/topspin3.6.3/exp/stan/nmr/lists/base\_info/...

[..]

Every non-Bruker file in /opt/topspin3.6.3/exp/stan/nmr/lists/vd/ will be saved:

now extracting from dir /opt/topspin3.6.3/exp/stan/nmr/lists/vd/...

Every non-Bruker file in /opt/topspin3.6.3/exp/stan/nmr/lists/vp/ will be saved:

now extracting from dir /opt/topspin3.6.3/exp/stan/nmr/lists/vp/...

Every non-Bruker file in /opt/topspin3.6.3/exp/stan/nmr/lists/vt/ will be saved:

now extracting from dir /opt/topspin3.6.3/exp/stan/nmr/lists/vt/...

Every non-Bruker file in /opt/topspin3.6.3/exp/stan/nmr/lists/wave/ will be saved:

now extracting from dir /opt/topspin3.6.3/exp/stan/nmr/lists/wave/...

Now writing zip file

/opt/topspin3.6.3/nmr\_backup/extractedDirectories.zip7...

Every non-Bruker file in /opt/topspin3.6.3/exp/stan/nmr/par/ will be saved:

```
now extracting from dir /opt/topspin3.6.3/exp/stan/nmr/par/...

Now writing zip file
/opt/topspin3.6.3/nmr_backup/extractedDirectories.zip8...

Every non-Bruker file in /opt/topspin3.6.3/exp/stan/nmr/py/ will be saved:
now extracting from dir /opt/topspin3.6.3/exp/stan/nmr/py/...

Now writing zip file
/opt/topspin3.6.3/nmr_backup/extractedDirectories.zip9...

Now writing zip file
/opt/topspin3.6.3/nmr_backup/nmr_backup_sermnadmin_20210812-1137.zip...

#####

Backing up of NMR data was successful!
The backup file

/opt/topspin3.6.3/nmr_backup/nmr_backup_sermnadmin_20210812-1137.zip

has been generated!

#####
```

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## exp/stan/nmr/lists/ (multiples carpetes)

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```
BIS file for magnet no 286 / cryostat 5051
```

Note:

manually creating a BIS file should be done for OLD systems only (e.g. no BIS available for download).

Do not attempt to use this function for systems with active cooling and/or liquefaction technology!

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```
BIS Parameter Definition
```

```
This form must be used for NMR magnets only.
```

```
Please do this very carefully.
```

```
Parameters should be entered based on magnet manual and type label. Entering wrong data might result in erroneous indications and/or false alarms in system monitoring.
```

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