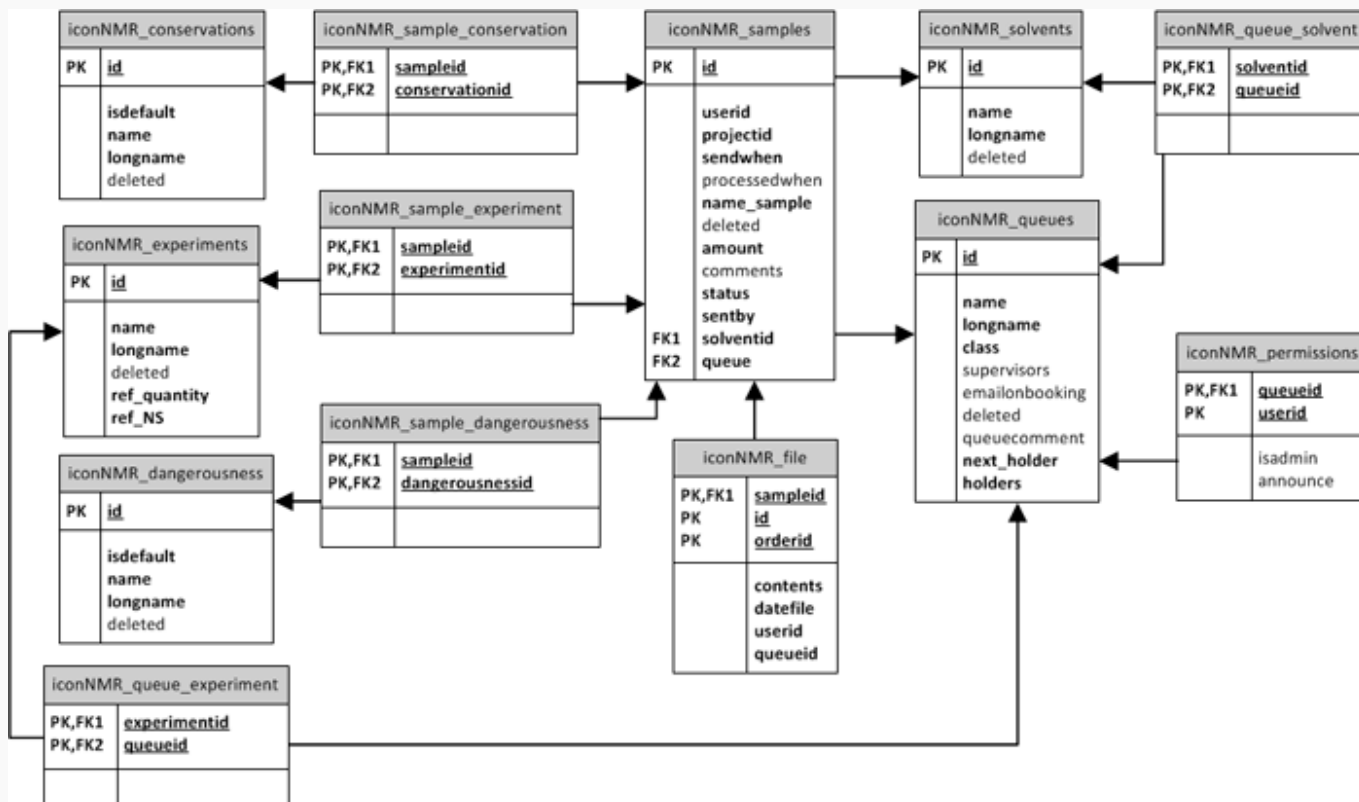
Cua de Rutina (Mòdul) - IconNMR

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```
# iconnmr external setup file
#
# dual experiment test file - default parameters modified
#
USER SeRMN
HOLDER 33
NAME ext-set-test-003
SOLVENT acetone
#
# begin of 1st experiment
#
EXPNO 1
NO_SUBMIT
PARAMETERS sw,20,olp,0,ns,512
EXPERIMENT teo_proto
TITLE ext-set-test-004\n 1D-1H-experiment
#
# end of 1st experiment
#
# begin of 2nd experiment
#
EXPNO 2
NO_SUBMIT
PARAMETERS ns,32k
EXPERIMENT teo_carboni
TITLE ext-set-test-00\n 1D-13C-experiment
#
# end of 2nd experiment
#
# end of holder definition
END
```

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```
# iconnmr external setup file
#
# three users

# ##### BEGIN USER
USER SeRMN

# ##### BEGIN SAMPLE (HOLDER)
```

```
HOLDER 31
DELETE
NAME user1-sample1
SOLVENT cdcl3
##### BEGIN EXPERIMENT
PARAMETERS ns,128,swp,15,td,8192
EXPERIMENT teo_proto
TITLE user1-sample1-exp1
##### END SAMPLE (HOLDER)
END
```

```
##### BEGIN USER
USER avirgili
```

```
##### BEGIN SAMPLE (HOLDER)
HOLDER 32
DELETE
NAME user2-sample1
SOLVENT dmso
##### BEGIN EXPERIMENT
PARAMETERS ns,128,swp,15,td,8192
EXPERIMENT teo_proto
TITLE user2-sample1-exp1
##### BEGIN EXPERIMENT
EXPERIMENT teo_carboni
TITLE user2-sample1-exp2
##### END SAMPLE (HOLDER)
END
```

```
##### BEGIN SAMPLE (HOLDER)
HOLDER 33
DELETE
NAME user2-sample2
SOLVENT dmso
##### BEGIN EXPERIMENT
PARAMETERS ns,128,swp,15,td,8192
EXPERIMENT teo_proto
TITLE user2-sample2-exp1
##### BEGIN EXPERIMENT
EXPERIMENT teo_carboni
TITLE user2-sample2-exp2
##### END SAMPLE (HOLDER)
END
```

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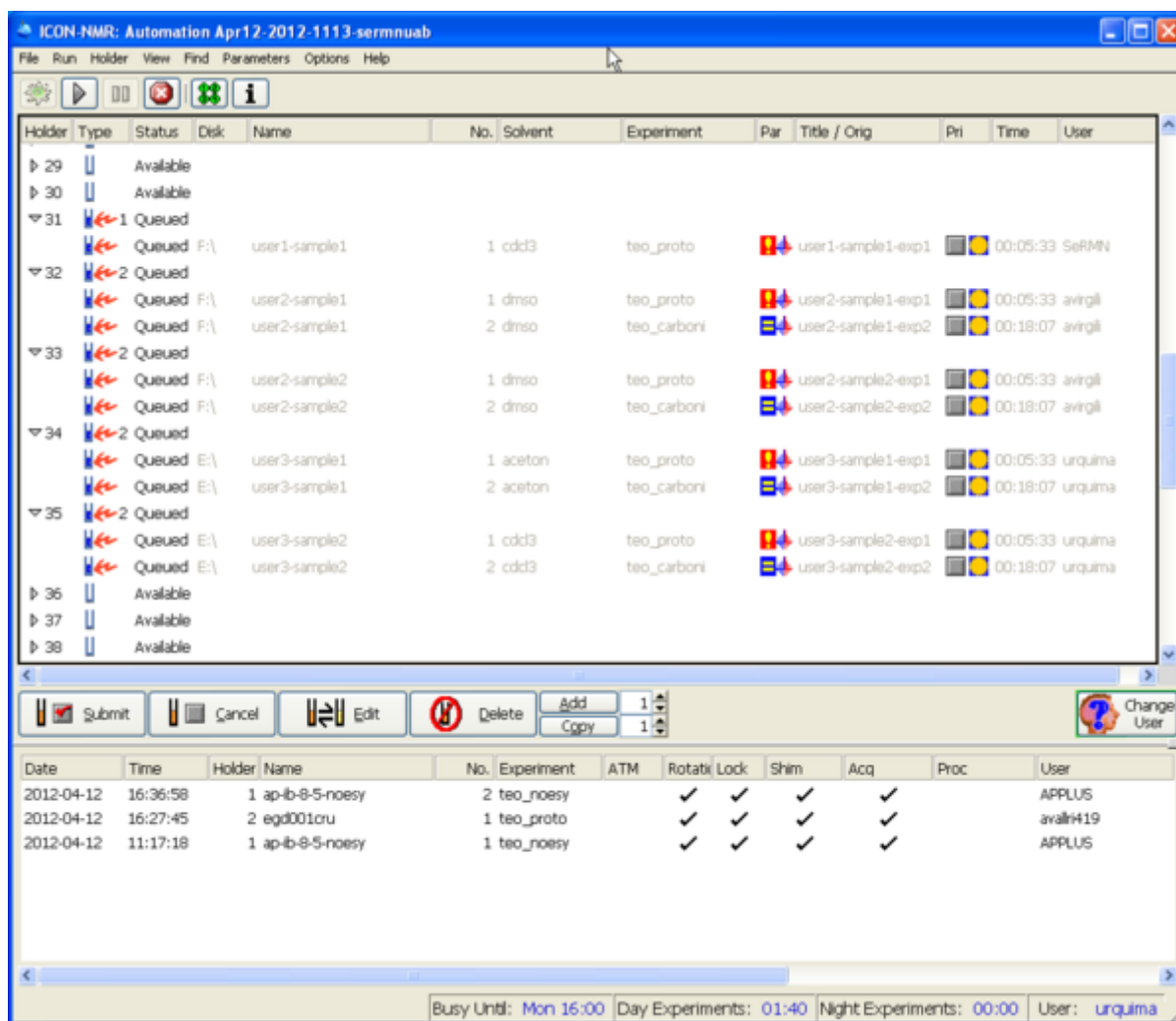
Change User

Busy Until: No Jobs! Day Experiments: 00:00 Night Experiments: 00:00 User: albert

TD 16384

NS 32

SW 10.9592 [ppm]

O1P 4.892

D1 1

OK

Index	State	Path	File Name	Size	Owner	Permissions	Mod Time	Mod User
▷ 29	Available							
▷ 30	Available							
▽ 31	1 Queued							
	Queued	D:\	ext-set-test-001	1 cdd3	teo_proto	ext-set-tes	00:02:21	sermn
▽ 32	1 Queued							
	Queued	D:\	ext-set-test-002	1 cdd3	teo_proto			sermn
▽ 33	3 Queued							
	Queued	D:\	ext-set-test-003	1 acetone	teo_proto	ext-set-tes 1D-1H-exp	00:02:21	sermn
	Available	D:\	ext-set-test-003	2 acetone	teo_carboni	ext-set-tes 1D-13C-ex		sermn
	Queued	D:\	ext-set-test-003	3 acetone	teo_proto		00:02:21	sermn
	Available	D:\	ext-set-test-003	4 acetone	teo_carboni	ext-set-tes 1D-1H-exp	02:01:01	sermn
▷ 34	Available							
▷ 35	Available							
▷ 36	Available							

Index	Status	Path	File Name	Size	Format	Thumbnail	Preview	Time	Owner
30	Available								
31	Available								
	Available	D:\ext-set-test-001	1 cdc3	teo_proto		ext-set-test-001			SeRMN
32	Available								
	Available	D:\ext-set-test-002	1 cdc3	teo_proto		ext-set-test-1H-experin		00:05:33	SeRMN
33	Available								
	Available	D:\ext-set-test-003	1 acetone	teo_proto		ext-set-test-1D-1H-exp			SeRMN
	Available	D:\ext-set-test-003	2 acetone	teo_carboni		ext-set-test-1D-13C-ex			SeRMN
	Available	D:\ext-set-test-003	3 acetone	teo_proto		ext-set-test-1D-1H-exp		00:22:59	SeRMN
	Available	D:\ext-set-test-003	4 acetone	teo_carboni		ext-set-test-1D-13C-ex		00:01:16	SeRMN
34	Available								
35	Available								

The screenshot displays the Bruker TopSpin software interface. The 'Acquisition Parameters' dialog box is open, showing the following parameters:

Parameter	Value	Unit
TD	8192	
NS	128	
SW	10.9592	[ppm]
O1P	4.892	
D1	1	

The 'OK' button is located at the bottom of the dialog box. The background shows the 'Acquisition Parameters' table with columns for 'Name', 'Value', 'Unit', and 'Comment'. The table lists parameters for 'ext-set-test-001' and 'ext-set-test-002'.

sermn_wiki:userpages:marta:cu_a_rutina https://sermn.uab.cat/wiki/doku.php?id=sermn_wiki:userpages:marta:cu_a_rutina&rev=1337856895

	Status	Path	Set Name	Count	Substance	Protocol	Acquisition Parameters	Time	Operator
	Available	D:\	ext-set-test-001	1	cdd3	teo_proto	ext-set-test-tes 1D-1H - cc		SerMN
	Available	D:\	ext-set-test-001	2	cdd3	teo_proto	ext-set-test-tes 1D-1H - cc		SerMN
	Available								
	Available	D:\	ext-set-test-002	1	cdd3	teo_proto	ext-set-test-tes 1D 1H - cc	00:05:33	SerMN
	Available								
	Available	D:\	ext-set-test-003	1	acetone	teo_proto	ext-set-test-tes 1D 1H - ac		SerMN
	Available	D:\	ext-set-test-003	2	acetone	teo_carboni	ext-set-test-tes 1D 13C - a		SerMN
	Available	D:\	ext-set-test-003	3	acetone	teo_proto	ext-set-test-tes 1D 1H - ac		SerMN
	Available	D:\	ext-set-test-003	4	acetone	teo_carboni	ext-set-test-tes 1D 13C - a		SerMN
	Available								
	Available	D:\	ext-set-test-004	1	acetone	teo_proto	<data - hor 1D 1H - ac	00:22:59	SerMN
	Available	D:\	ext-set-test-004	2	acetone	teo_carboni	<data - hor 1D-13C - a	00:01:16	SerMN

From: <https://sermn.uab.cat/wiki/> - **SeRMN Wiki**

Permanent link:
https://sermn.uab.cat/wiki/doku.php?id=sermn_wiki:userpages:marta:cu_a_rutina&rev=1337856895

Last update: **2012/05/24 12:54**