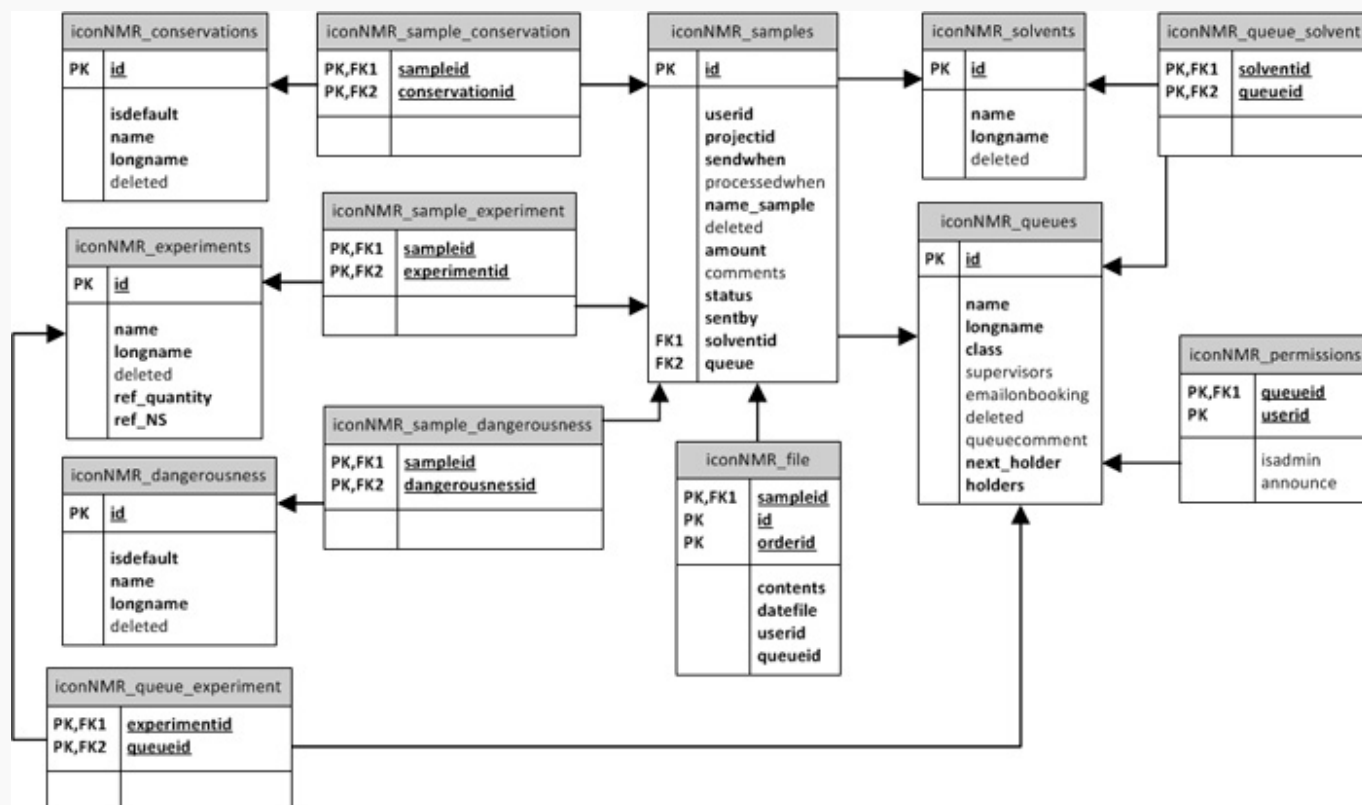
Cua de Rutina (Mòdul) - IconNMR

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```
# iconnmr external setup file
#
# dual experiment test file - default parameters modified
#
USER SeRMN
HOLDER 33
NAME ext-set-test-003
SOLVENT acetone
#
# begin of 1st experiment
#
EXPNO 1
NO_SUBMIT
PARAMETERS sw,20,olp,0,ns,512
EXPERIMENT teo_proto
TITLE ext-set-test-004\n 1D-1H-experiment
#
# end of 1st experiment
#
# begin of 2nd experiment
#
EXPNO 2
NO_SUBMIT
PARAMETERS ns,32k
EXPERIMENT teo_carboni
TITLE ext-set-test-00\n 1D-13C-experiment
#
# end of 2nd experiment
#
# end of holder definition
END
```

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```
# iconnmr external setup file
#
# three users
```

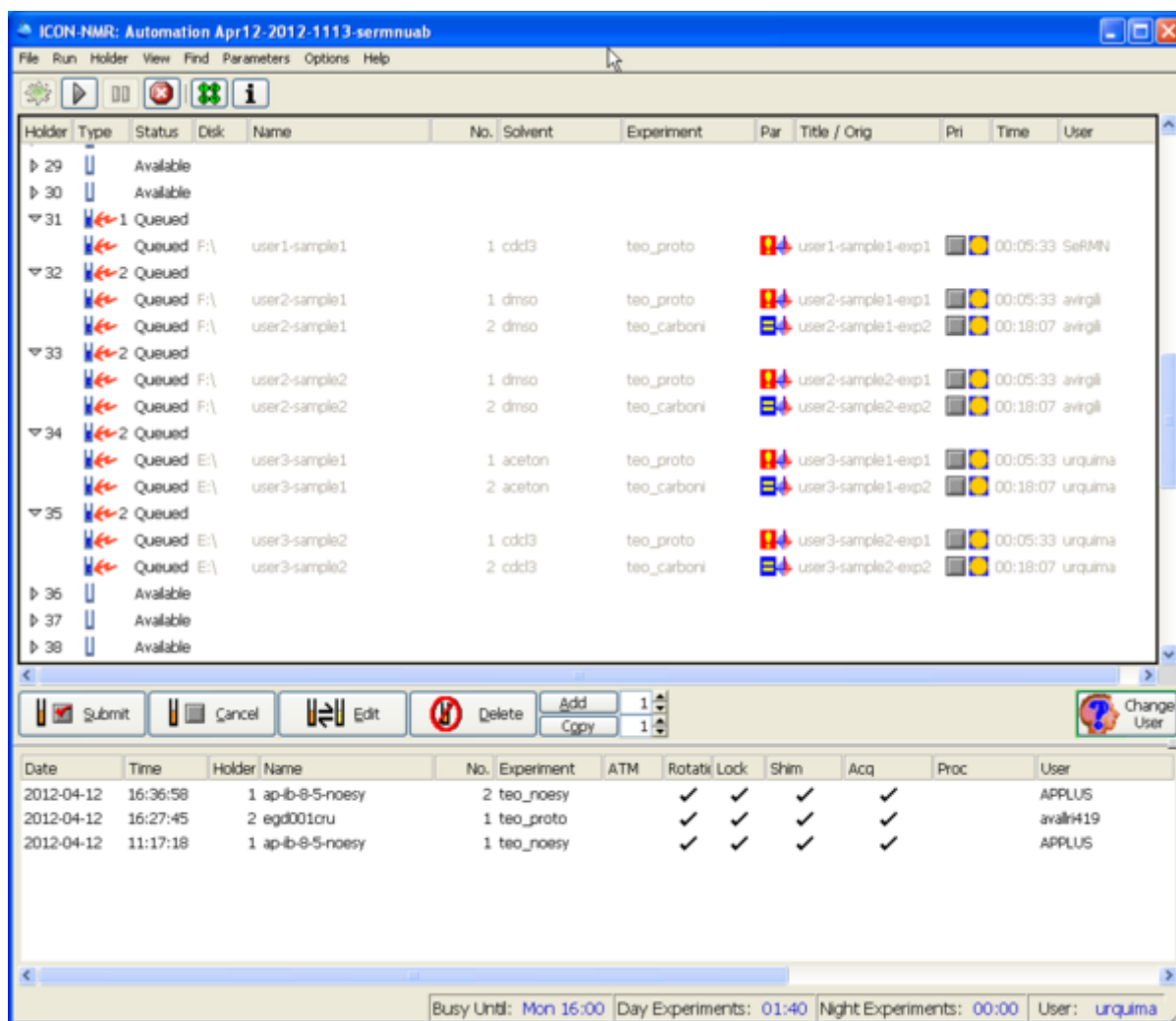
```
# ##### BEGIN USER
USER SeRMN

# ##### BEGIN SAMPLE (HOLDER)
HOLDER 31
DELETE
NAME user1-sample1
SOLVENT cdcl3
# ##### BEGIN EXPERIMENT
PARAMETERS ns,128,swp,15,td,8192
EXPERIMENT teo_proto
TITLE user1-sample1-exp1
##### END SAMPLE (HOLDER)
END

##### BEGIN USER
USER avirgili

##### BEGIN SAMPLE (HOLDER)
HOLDER 32
DELETE
NAME user2-sample1
SOLVENT dmso
##### BEGIN EXPERIMENT
PARAMETERS ns,128,swp,15,td,8192
EXPERIMENT teo_proto
TITLE user2-sample1-exp1
##### BEGIN EXPERIMENT
EXPERIMENT teo_carboni
TITLE user2-sample1-exp2
##### END SAMPLE (HOLDER)
END

##### BEGIN SAMPLE (HOLDER)
HOLDER 33
DELETE
NAME user2-sample2
SOLVENT dmso
##### BEGIN EXPERIMENT
PARAMETERS ns,128,swp,15,td,8192
EXPERIMENT teo_proto
TITLE user2-sample2-exp1
##### BEGIN EXPERIMENT
EXPERIMENT teo_carboni
TITLE user2-sample2-exp2
##### END SAMPLE (HOLDER)
END

##### BEGIN USER
```

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Change User

Busy Until: No Jobs! Day Experiments: 00:00 Night Experiments: 00:00 User: albert

TD 16384

NS 32

SW 10.9592 [ppm]

O1P 4.892

D1 1

OK

Task ID	Status	Path	Test Name	Count	Substrate	Protocol	Test Name	Time	Operator
29	Available								
30	Available								
31	1 Queued								
	Queued	D:\	ext-set-test-001	1	cdd3	teo_proto	ext-set-tes	00:02:21	sermn
32	1 Queued								
	Queued	D:\	ext-set-test-002	1	cdd3	teo_proto			sermn
33	3 Queued								
	Queued	D:\	ext-set-test-003	1	acetone	teo_proto	ext-set-tes 1D-1H-exp	00:02:21	sermn
	Available	D:\	ext-set-test-003	2	acetone	teo_carboni	ext-set-tes 1D-13C-ex		sermn
	Queued	D:\	ext-set-test-003	3	acetone	teo_proto		00:02:21	sermn
	Available	D:\	ext-set-test-003	4	acetone	teo_carboni	ext-set-tes 1D-1H-exp	02:01:01	sermn
34	Available								
35	Available								
36	Available								

Index	Status	Path	Size	File Type	Protocol	Thumbnail	File Name	Duration	Source
30	Available								
31	Available								
	Available	D:\ ext-set-test-001	1 cdd3	teo_proto		ext-set-tes			SeRMN
32	Available								
	Available	D:\ ext-set-test-002	1 cdd3	teo_proto		ext-set-tes 1H-experin		00:05:33	SeRMN
33	Available								
	Available	D:\ ext-set-test-003	1 aceton	teo_proto		ext-set-tes 1D-1H-exp			SeRMN
	Available	D:\ ext-set-test-003	2 aceton	teo_carboni		ext-set-tes 1D-13C-ex			SeRMN
	Available	D:\ ext-set-test-003	3 aceton	teo_proto		ext-set-tes 1D-1H-exp		00:22:59	SeRMN
	Available	D:\ ext-set-test-003	4 aceton	teo_carboni		ext-set-tes 1D-13C-ex		00:01:16	SeRMN
34	Available								
35	Available								

The screenshot shows the Bruker TopSpin software interface. The 'Acquisition Parameters' dialog box is open, displaying the following parameters:

Parameter	Value	Unit	Comment
TD	8192		
NS	128		
SW	10.9592	[ppm]	
O1P	4.892		
D1	1		

The 'OK' button is visible at the bottom of the dialog box. The background shows the 'Acquisition Parameters' table with columns for 'Name', 'Value', 'Unit', and 'Comment'.

sermn_wiki:userpages:marta:cu_a_rutina https://sermn.uab.cat/wiki/doku.php?id=sermn_wiki:userpages:marta:cu_a_rutina&rev=1337857476

	Available	D:\	ext-set-test-001	1	cdd3	teo_proto	 ext-set-test-1D-1H - cc 	SeRMN
	Available	D:\	ext-set-test-001	2	cdd3	teo_proto	 ext-set-test-1D-1H - cc 	SeRMN
	Available							
	Available	D:\	ext-set-test-002	1	cdd3	teo_proto	 ext-set-test-1D 1H - cc  00:05:33	SeRMN
	Available							
	Available	D:\	ext-set-test-003	1	acetone	teo_proto	 ext-set-test-1D 1H - ac 	SeRMN
	Available	D:\	ext-set-test-003	2	acetone	teo_carboni	 ext-set-test-1D 13C - a 	SeRMN
	Available	D:\	ext-set-test-003	3	acetone	teo_proto	 ext-set-test-1D 1H - ac 	SeRMN
	Available	D:\	ext-set-test-003	4	acetone	teo_carboni	 ext-set-test-1D 13C - a 	SeRMN
	Available							
	Available	D:\	ext-set-test-004	1	acetone	teo_proto	 <data - hor 1D 1H - ac  00:22:59	SeRMN
	Available	D:\	ext-set-test-004	2	acetone	teo_carboni	 <data - hor 1D-13C - a  00:01:16	SeRMN

From: <https://sermn.uab.cat/wiki/> - **SeRMN Wiki**

Permanent link:
https://sermn.uab.cat/wiki/doku.php?id=sermn_wiki:userpages:marta:cu_a_rutina&rev=1337857476

Last update: **2012/05/24 13:04**