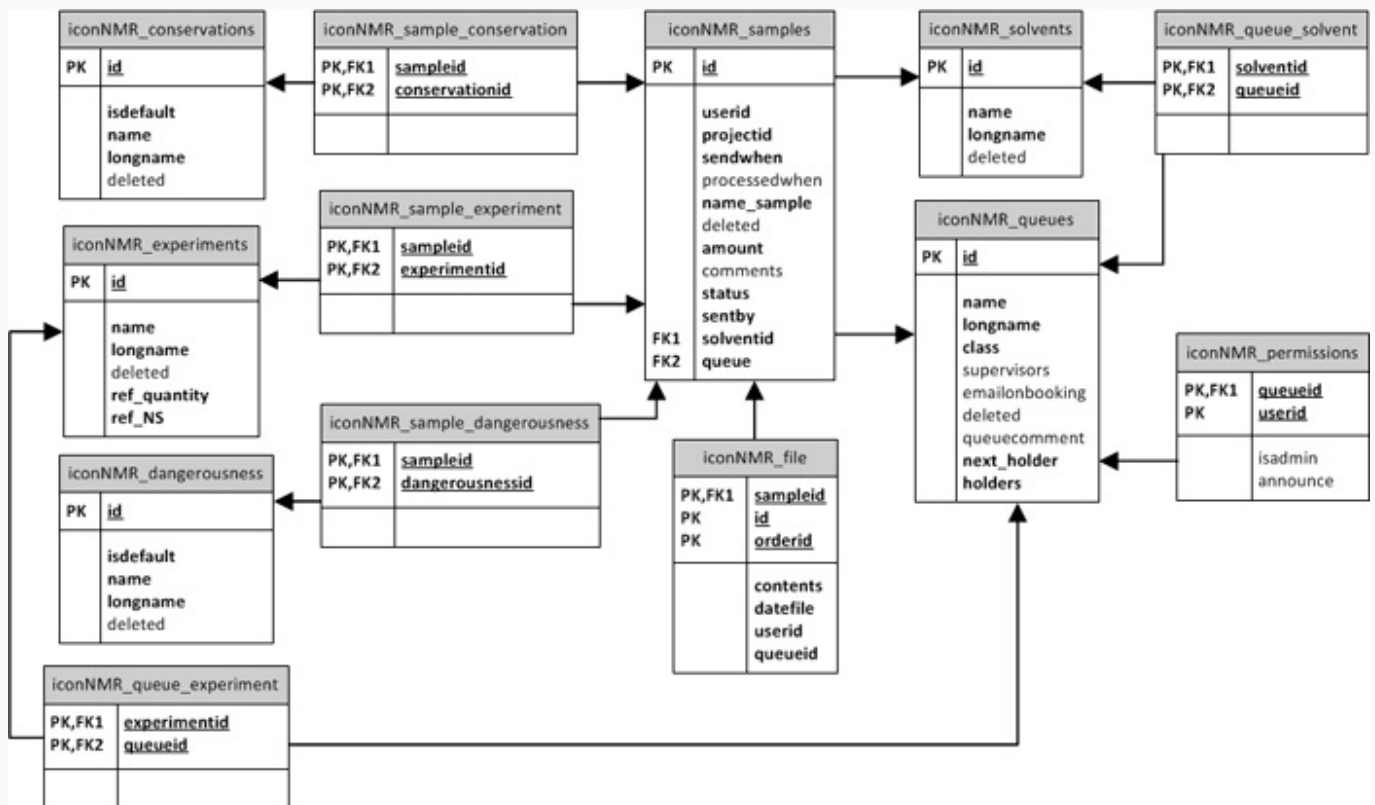
# Cua de Rutina (Mòdul) - IconNMR

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```
# iconnmr external setup file
#
# dual experiment test file - default parameters modified
#
USER SeRMN
HOLDER 33
NAME ext-set-test-003
SOLVENT acetone
#
# begin of 1st experiment
#
EXPNO 1
NO_SUBMIT
PARAMETERS sw,20,olp,0,ns,512
EXPERIMENT teo_proto
TITLE ext-set-test-004\n 1D-1H-experiment
#
# end of 1st experiment
#
# begin of 2nd experiment
#
EXPNO 2
NO_SUBMIT
PARAMETERS ns,32k
EXPERIMENT teo_carboni
TITLE ext-set-test-00\n 1D-13C-experiment
#
# end of 2nd experiment
#
# end of holder definition
END
```

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```
# iconnmr external setup file
#
# three users
```

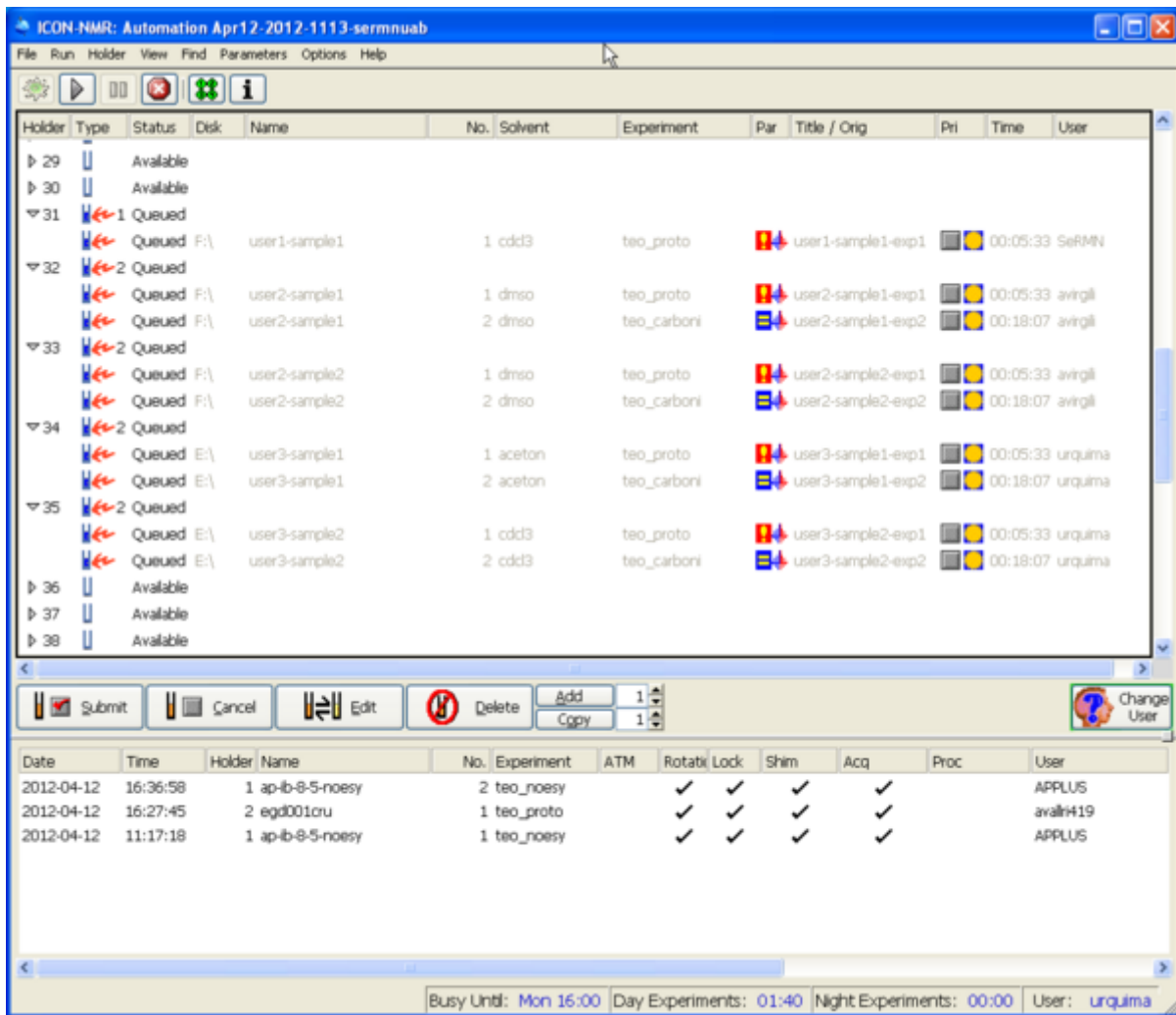
```
# #### BEGIN USER
USER SeRMN

# #### #### BEGIN SAMPLE (HOLDER)
HOLDER 31
DELETE
NAME user1-sample1
SOLVENT cdcl3
# #### #### #### BEGIN EXPERIMENT
PARAMETERS ns,128,swp,15,td,8192
EXPERIMENT teo_proto
TITLE user1-sample1-exp1
#### #### END SAMPLE (HOLDER)
END

#### BEGIN USER
USER avirgili

#### #### BEGIN SAMPLE (HOLDER)
HOLDER 32
DELETE
NAME user2-sample1
SOLVENT dmso
#### #### #### BEGIN EXPERIMENT
PARAMETERS ns,128,swp,15,td,8192
EXPERIMENT teo_proto
TITLE user2-sample1-exp1
#### #### #### BEGIN EXPERIMENT
EXPERIMENT teo_carboni
TITLE user2-sample1-exp2
#### #### END SAMPLE (HOLDER)
END

#### #### BEGIN SAMPLE (HOLDER)
HOLDER 33
DELETE
NAME user2-sample2
SOLVENT dmso
#### #### #### BEGIN EXPERIMENT
PARAMETERS ns,128,swp,15,td,8192
EXPERIMENT teo_proto
TITLE user2-sample2-exp1
#### #### #### BEGIN EXPERIMENT
EXPERIMENT teo_carboni
TITLE user2-sample2-exp2
#### #### END SAMPLE (HOLDER)
END

#### BEGIN USER
```

```
USER urquima

#### #### BEGIN SAMPLE (HOLDER)
HOLDER 34
DELETE
NAME user3-sample1
SOLVENT acetone
#### #### #### BEGIN EXPERIMENT
PARAMETERS ns,128,swp,15,td,8192
EXPERIMENT teo_proto
TITLE user3-sample1-exp1
#### #### #### BEGIN EXPERIMENT
EXPERIMENT teo_carboni
TITLE user3-sample1-exp2
#### #### END SAMPLE (HOLDER)
END

#### #### BEGIN SAMPLE (HOLDER)
HOLDER 35
DELETE
NAME user3-sample2
SOLVENT cdcl3
#### #### #### BEGIN EXPERIMENT
PARAMETERS ns,128,swp,15,td,8192
EXPERIMENT teo_proto
TITLE user3-sample2-exp1
#### #### #### BEGIN EXPERIMENT
EXPERIMENT teo_carboni
TITLE user3-sample2-exp2
#### #### END SAMPLE (HOLDER)
END
```

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The screenshot shows the Bruker TopSpin software interface. The main window displays a list of experiments with columns for Holder, Type, Status, Disk, Name, No., Solvent, Experiment, Par, Title, Orig. Prt, Time, and User. The experiments listed are mostly 'Finished' and include details like '1 CDC3', 'teo\_proto', and 'Synthesis'. A status window titled 'Feb01-2012-1223- Albert Status' is overlaid on the right, showing a grid of colored circles representing the status of various channels or parameters. Below the main window is a toolbar with buttons for Submit, Cancel, Edit, Delete, Add, and Copy. At the bottom, there is a detailed log table with columns for Date, Time, Holder, Name, No., Experiment, ATM, Rotafix, Lock, Shim, Acq, Proc, User, Disk, and TR Remarks. The status bar at the bottom indicates 'Busy Until: No Jobs!', 'Day Experiments: 00:00', 'Night Experiments: 00:00', and 'User: Albert'.

A dialog box with a blue icon in the top left corner. It contains five input fields with the following labels and values: TD (16384), NS (32), SW (10.9592 [ppm]), O1P (4.892), and D1 (1). An 'OK' button is located at the bottom left, and a mouse cursor is pointing at the right side of the dialog.

▷ 29		Available										
▷ 30		Available										
▽ 31		1 Queued										
		Queued	D:\	ext-set-test-001	1	cdd3	teo_proto		ext-set-tes		00:02:21	sermn
▽ 32		1 Queued										
		Queued	D:\	ext-set-test-002	1	cdd3	teo_proto					sermn
▽ 33		3 Queued										
		Queued	D:\	ext-set-test-003	1	acetone	teo_proto		ext-set-tes		00:02:21	sermn
		Available	D:\	ext-set-test-003	2	acetone	teo_carboni		ext-set-tes			sermn
		Queued	D:\	ext-set-test-003	3	acetone	teo_proto				00:02:21	sermn
		Available	D:\	ext-set-test-003	4	acetone	teo_carboni		ext-set-tes		02:01:01	sermn
▷ 34		Available										
▷ 35		Available										
▷ 36		Available										
▷ 37		Available										

▷ 30		Available										
▽ 31		Available										
		Available	D:\	ext-set-test-001	1	cdd3	teo_proto		ext-set-tes			SeRMN
▽ 32		Available										
		Available	D:\	ext-set-test-002	1	cdd3	teo_proto		ext-set-tes		00:05:33	SeRMN
▽ 33		Available										
		Available	D:\	ext-set-test-003	1	aceton	teo_proto		ext-set-tes			SeRMN
		Available	D:\	ext-set-test-003	2	aceton	teo_carboni		ext-set-tes			SeRMN
		Available	D:\	ext-set-test-003	3	aceton	teo_proto		ext-set-tes		00:22:59	SeRMN
		Available	D:\	ext-set-test-003	4	aceton	teo_carboni		ext-set-tes		00:01:16	SeRMN
▷ 34		Available										
▷ 35		Available										

▷ 30		Available										
▽ 31		Available										
		Available	D:\	ext-set-test-001	1	cdd3	teo_proto		ext-set			SeRMN
▽ 32		Available										
		Available	D:\	ext-set-test-002	1	cdd3	teo_proto		ext-set		00:05:33	SeRMN
▽ 33		Available										
		Available	D:\	ext-set-test-003	1	aceton	teo_proto					
		Available	D:\	ext-set-test-003	2	aceton	teo_carboni					
		Available	D:\	ext-set-test-003	3	aceton	teo_proto					
		Available	D:\	ext-set-test-003	4	aceton	teo_carboni					
▷ 34		Available										
▷ 35		Available										

TD	<input type="text" value="8192"/>	
NS	<input type="text" value="128"/>	
SW	<input type="text" value="10.9592"/>	[ppm]
O1P	<input type="text" value="4.892"/>	
D1	<input type="text" value="1"/>	
<input type="button" value="OK"/>		

▽ 31	Available							
	Available	D:\ ext-set-test-001	1	cdd3	teo_proto		ext-set-tes 1D-1H - cc	SeRMN
	Available	D:\ ext-set-test-001	2	cdd3	teo_proto		ext-set-tes 1D-1H - cc	SeRMN
▽ 32	Available							
	Available	D:\ ext-set-test-002	1	cdd3	teo_proto		ext-set-tes 1D 1H - cc	00:05:33 SeRMN
▽ 33	Available							
	Available	D:\ ext-set-test-003	1	aceton	teo_proto		ext-set-tes 1D 1H - ac	SeRMN
	Available	D:\ ext-set-test-003	2	aceton	teo_carboni		ext-set-tes 1D 13C - a	SeRMN
	Available	D:\ ext-set-test-003	3	aceton	teo_proto		ext-set-tes 1D 1H - ac	SeRMN
	Available	D:\ ext-set-test-003	4	aceton	teo_carboni		ext-set-tes 1D 13C - a	SeRMN
▽ 34	Available							
	Available	D:\ ext-set-test-004	1	aceton	teo_proto		<data - hor 1D 1H - ac	00:22:59 SeRMN
	Available	D:\ ext-set-test-004	2	aceton	teo_carboni		<data - hor 1D-13C - a	00:01:16 SeRMN
▷ 35	Available							
▷ 36	Available							

From: <https://sermn.uab.cat/wiki/> - **SeRMN Wiki**

Permanent link: [https://sermn.uab.cat/wiki/doku.php?id=sermn\\_wiki:userpages:marta:cua\\_rutina&rev=1338454396](https://sermn.uab.cat/wiki/doku.php?id=sermn_wiki:userpages:marta:cua_rutina&rev=1338454396)

Last update: **2012/05/31 10:53**