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The screenshot shows the 'SeRMN Instrument Booking Tool UAB' interface. At the top, it says 'You are logged as: usuari.normal'. The main heading is 'Select an instrument to manage your bookings'. Below this is a table with columns 'Instrument', 'Description', and 'Location'. The table lists three instruments: '250 MHz - AUTOSERVEI', '400 MHz', and 'Dynamic Nuclear Polarizer'. To the left is a navigation menu with options like 'Main Menu', 'Instruments & Queues', 'My bookings', 'My account', 'Change password', 'Logout', 'IconNMR module', 'My samples', 'Help contents', 'About', and 'Help'. Below the instrument table is another section 'Select a queue to send a sample - IconNMR' with a table showing 'Cua_1' and its description.

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System managed by **Bumblebee** version 1.3.3. This installation of Bumblebee currently manages 294 users, 57 projects, 17 instruments and 9543 bookings. Running under Linux (2.6.32-5-686), Apache (2.2.16), PHP (5.3.3-7squeeze3) and MySQL (5.1.61-0+squeeze1).

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Solvent	AcCapil (Acetone-d6 in Capillary Tube)
Sample name	
Amount (mg)	50

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- [00-UserQueue](#)

[Return to main menu](#)

System managed by **Bumblebee** version 1.3.1, released under the GNU GPL.
 This installation of Bumblebee currently manages 294 users, 57 projects, 17 instruments and 9543 bookings.
 Running under Linux (2.6.32-5-686), Apache (2.2.16), PHP (5.3.3-1ubuntu5) and MySQL (5.1.40-1ubuntu10).
 Email the **system administrator** for help.
Booking information Copyright © 2012 Servei de Resonància Magnètica Nuclear, UAB

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Sample could not be made:
Some values entered into the form are not valid and should be highlighted in the form below. Please check your data entry and try again.

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- **Test 1** some important information about the queue can be here.
- **Test 2** some important information about the queue can be here.

User	296 (UsuariNormal)
Recorded by	296 (UsuariNormal)
Queue	1 (Cua_1)
Sent Date	2012-05-24 10:50
My Projects	Teo - SeRMN (T063300 / PK122933 - 2010)
Solvent	AcCapil (Acetone-d6 in Capillary Tube)
Sample name	<input type="text"/>
Amount (mg)	50

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Experiments and properties

Experiments (Press CTRL to multiple choice)

(Select one or more options)
Teo Carboni
Teo H₂O
Teo Noesy

Dangerousness (Press CTRL to multiple choice)

(Select one or more options)
(Unknown)
Medium toxic level
Inflammable description

Conservation (Press CTRL to multiple choice)

(Select one or more options)
Temperature between 20-30°C
Freeze -20°C

Send sample

Queue supervisors

- [00-UserQueue](#)

[Return to main menu](#)

System managed by **Bumblebee** version 1.3.1, released under the GNU GPL.
This installation of Bumblebee currently manages 294 users, 57 projects, 17 instruments and 9543 bookings.
Running under Linux (2.6.32-5-686), Apache (2.2.16), PHP (5.3.3-1ubuntu5) and MySQL (5.1.40-1ubuntu1).
Email the **system administrator** for help.

Booking information Copyright © 2012 Servei de Resonància Magnètica Nuclear, UAB

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- **Test** some important information about the queue can be here.
- **Test 2** some important information about the queue can be here.

User 296 (UsuariNormal)

Recorded by 296 (UsuariNormal)

Created by 296 (UsuariNormal)

Sent Date 2012-05-24 10:52

My Projects Teo - SeRMN (T063300 / PK122933 - 2010)

Solvent AcCapil (Acetone-d6 in Capillary Tube)

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The screenshot shows a web form for booking samples. It includes several dropdown menus and a 'Send sample' button. The interface is annotated with yellow callout boxes containing instructions in Catalan. The callouts are: 1. A blue box pointing to a dropdown menu with the text 'Els camps a emplenar apareixen en vermell perquè els puguis identificar ràpidament.' (The fields to be filled appear in red so you can identify them quickly). 2. A red box pointing to a dropdown menu with the text 'Els valors seleccionats apareixen amb un color diferent perquè els puguis distingir clarament.' (The selected values appear with a different color so you can distinguish them clearly). 3. A blue box pointing to the 'Send sample' button with the text 'Quan hagi corregit els errors, prem el botó [Send sample] per registrar la teva petició.' (When you have corrected the errors, press the [Send sample] button to register your request). 4. A blue box pointing to a link at the bottom with the text 'Però si canvies d'opinió i no vols sol·licitar res, clica l'enllaç "Return to main menu" per tornar a la pantalla principal.' (But if you change your mind and do not want to request anything, click the link "Return to main menu" to return to the main screen).

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System managed by **Bumblebee** version 1.3.1, released under the GNU GPL.
This installation of Bumblebee currently manages 294 users, 57 projects, 11 instruments and 9543 bookings.
Running under Linux (2.6.32-5-686), Apache (2.2.16), PHP (5.3.3) and MySQL (5.1.40).
Email the **system administrator** for help.
Booking information Copyright © 2012 Servei de Resonància Magnètica Nuclear, UAB

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English

Main Menu

- Instruments & Queues
- My bookings
- My account
- Change password
- Logout
- IconNMR module
- My samples
- Help contents
- About
- Help

Sample made

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- **Test** some important information about the queue can be here.
- **Test 2** some important information about the queue can be here.

Sample ID	13
User	296 (UsuariNormal)
Recorded by	296 (UsuariNormal)
Queue	1 (Cua_1)
Sent Date	2012-05-24 10:52
My Projects	Teo - SeRMN (T063300 / PK122933 - 2010)

Solvent	AcCapil (Acetone-d6 in Capillary Tube)
Sample name	sample for nmr
Amount (mg)	50

Comments

Experiments and properties

Experiment (Press CTRL to multiple choice)	(Select one or more options) Teo Carboni
--	---

Experiments and properties

Experiment (Press CTRL to multiple choice)	(Select one or more options) Teo Carboni Teo HMBC Teo Carb
--	---

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Dangerousness (Press CTRL to multiple choice)	(Select one or more options) (Unknown) Medium toxic level Inflammable description
---	--

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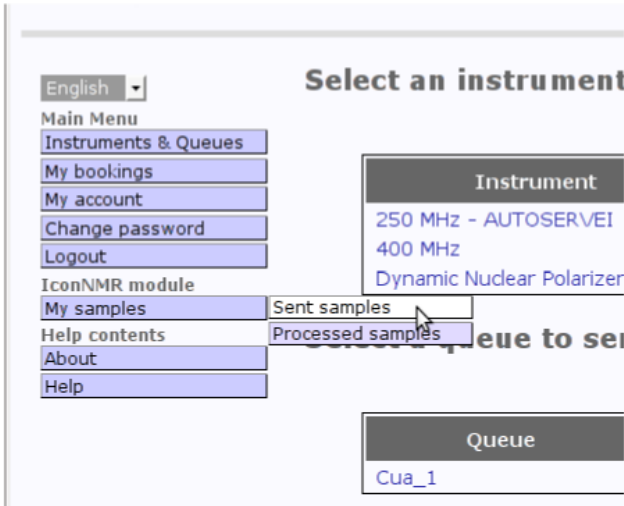
System managed by **Bumblebee** version 1.3.1, released under the GNU GPL. This installation of Bumblebee currently manages 294 users, 57 projects, 17 instruments and 9543 bookings. Running under Linux (2.6.32-5-686), Apache (2.2.16), PHP (5.3.3-1ubuntu3) with MySQL (5.1.40-1ubuntu3) and InfluxDB (0.9.0-1ubuntu1).

Email the **system administrator** for help.

Booking information Copyright © 2012 Servei de Resonància Magnètica Nuclear, UAB

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System managed by **Bumblebee** version 1.3.1, released under the GPL. This installation of Bumblebee currently manages 294 users, 57 projects, 17 instruments, 9543 bookings, sample submission, folios and requests. Running under Linux (2.6.32-5-686), Apache (2.2.16), PHP (5.3.3-7+squeezed-apache2handler-mode) with MySQL (5.1.61-0+squeezed1).

Email the **system administrator** for help.

Booking information Copyright © 2012 Servei de Resonància Magnètica Nuclear, UAB

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System managed by **Bumblebee** version 1.3.1, released under the [GNU GPL](#).
This installation of Bumblebee currently manages 294 users, 57 projects, 1744 samples, 1244 bookings, 3440 sample acquisition forms and requests.
Running under Linux (2.6.32-5-686), Apache (2.2.16), PHP (5.3.3-7+squeezeb, apache2handler mode) with MySQL (5.1.61-0+squeezel).

Email the **system administrator** for help.

Booking information Copyright © 2012 Servei de Resonància Magnètica Nuclear, UAB

SeRMN Instrument Booking Tool UAB

You are logged as: usuari.normal

English ▾

Main Menu

- Instruments & Queues
- My bookings
- My account
- Change password
- Logout
- IconNMR module
- My samples
- Help contents
- About
- Help

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- **Test** some important information about the queue can be here.
- **Test 2** some important information about the queue can be here.

Sample ID	13
User	296 (UsuariNormal)
Recorded by	296 (UsuariNormal)
Queue	1 (Cua_1)
Sent Date	2012-05-24 10:52
My Projects	Teo - SeRMN (T063300 / PK122933 - 2010)

Solvent	AcCapil (Acetone d6 in Capillary Tube)
Sample name	sample for nmr
Amount (mg)	50

Comments	
----------	--

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Experiments and properties

Experiment (Press CTRL to multiple choice)	(Select one or more options) Teo Carboni Teo HMBC Teo Noesy
Dangerousness (Press CTRL to multiple choice)	(Select one or more options) (Unknown) Medium toxic level Inflammable description
Conservation (Press CTRL to multiple choice)	(Select one or more options) Temperature between 20-30°C Freeze -20°C

Queue supervisors

- [00-UserQueue](#)

[Go to my samples](#)

System managed by **Bumblebee** version 1.3.1, released under the **GNU GPL**.
This installation of Bumblebee currently manages 294 users, 57 projects, 17 instruments, 9543 bookings, 2 sample submission forms and 14 requests.
Running under Linux (2.6.32-5-686), Apache (2.2.16), PHP (5.3.3-7+squeeze8, apache2handler mode) with MySQL (5.1.61-0+squeeze1).

Email the **system administrator** for help.

Booking information Copyright © 2012 Servei de Resonància Magnètica Nuclear, UAB

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Select an instrument to manage your bookings

- English
- Main Menu
 - Instruments & Queues
 - My bookings
 - My account
 - Change password
 - Logout
- IconNMR module
 - My samples
 - Help contents
 - About
 - Help

Instrument	Description	Location
250 MHz - AUTOSERVEI	Bruker Avance DRX-250 autoservei	SeRMN UAB
400 MHz	Bruker Avance III 400	SeRMN-UAB
Dynamic Nuclear Polarizer	Oxford Hypersense DNP	SeRMN UAB

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System managed by Bumblebee version 1.3.1, released under the GNU GPL... This installation of Bumblebee currently manages 294 users, 57 projects, 17 instruments, 9543 bookings, 2 sample submission forms and 19 requests.

Email the system administrator for help.

Booking information Copyright © 2012 Servei de Resonància Magnètica Nuclear, UAB

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- **Test** some important information about the queue can be here.
- **Test 2** some important information about the queue can be here.

Sample ID	13
User	296 (UsuariNormal)
Recorded by	296 (UsuariNormal)
Queue	1 (Cua_1)
Sent Date	2012-05-24 10:52
My Projects	Teo - SeRMN (T063300 / PK122933 - 2010)
Solvent	AcCapil (Acetone-d6 in Capillary Tube)
Sample name	sample for nmr
Amount (mg)	50
Comments	

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Update sample Delete sample Sample as new

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- **Test** some important information about the queue can be here.
- **Test 2** some important information about the queue can be here.

Sample ID: 13
 User: 296 (UsuariNormal)
 Recorded by: 296 (UsuariNormal)
 Queue: 1 (Cua_1)
 Sent Date: 2012-05-24 10:52
 My Projects: Teo - SeRMN (T063300 / PK122933 - 2010)

Are you sure to delete this sample? (this action cannot be undone)

Cancel OK

Experiments and properties

Experiment (Press CTRL to multiple choice): (Select one or more options)
 Teo Carboni
 Teo Noesy

Dangerousness (Press CTRL to multiple choice): (Select one or more options)
 Inflammable description

Conservation (Press CTRL to multiple choice): (Select one or more options)
 temperature between 20-30°C
 Freeze -20°C

Update sample Delete sample Sample as new

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You are logged in as UsuariNormal

English ▾

- Main Menu
- Instruments & Queues**
- My bookings
- My account
- Change password
- Logout
- IconNMR module
- My samples
- Help contents
- About
- Help

My sent samples (you can edit)

2012-05-18 16:07 to 2012-05-25 16:07 [< earlier](#) | [now](#) | [later >](#)

Sent Date	Sample name	Queue	Project	Sent by
2012-05-24 11:09	another sample for n	Cua_1	Teo - SeRMN	UsuariNormal
2012-05-23 11:40	Normal_1	Cua_1	Bioiberica	UsuariNormal

System managed by **Bumblebee** version 1.3.1, released under the **GNU GPL**.
 This installation of Bumblebee currently manages 294 users, 57 projects, 17 instruments, 9543 bookings, 2 sample submission forms and 19 requests.
 Running under Linux (2.6.32-5-686), Apache (2.2.16), PHP (5.3.3-7+squeezee8, apache2handler mode) with MySQL (5.1.61-0+squeezee1).

Email the **system administrator** for help.

Booking information Copyright © 2012 Servei de Resonància Magnètica Nuclear, UAB

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English ▾

Main Menu

- Instruments & Queues
- My bookings
- My account
- Change password
- Logout

IconNMR module

- My samples
- Processed samples
- Help contents
- About
- Help

Select an instrument to manage your bookings

Instrument	Description	Location
250 MHz - AUTOSERVEI	Bruker Avance DRX-250 autoservei	SeRMN UAB
400 MHz	Bruker Avance III 400	SeRMN-UAB
Dynamic Nuclear Polarizer	Oxford Hypersense DNP	SeRMN UAB

Queue to send a sample - IconNMR

Queue	Description
Cua_1	

System managed by **Bumblebee** version 1.3.1, released under the GPL license.
This installation of Bumblebee currently manages 294 users, 57 projects, 17 instruments, 9543 bookings, 2 sample submission forms and 19 requests.
Running under Linux (2.6.32-5-686), Apache (2.2.16), PHP (5.3.3-7+squeeze8, apache2handler mode) with MySQL (5.1.61-0+squeeze1).

Email the **system administrator** for help.

Booking information Copyright © 2012 Servei de Resonància Magnètica Nuclear, UAB

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English ▾

Main Menu

- Instruments & Queues
- My bookings
- My account
- Change password
- Logout

IconNMR module

- My samples
- Help contents
- About
- Help

My processed samples

2012-05-18 16:10 to 2012-05-25 16:10 [« earlier](#) | [now](#) | [later »](#)

Sent Date	Processed Date	Sample name	Queue	Project	Sent by
2012-05-23 11:40	2012-05-25 14:55	Normal_2	Cua_1	Bioiberica	UsuariNormal
2012-05-23 18:16	2012-05-25 14:55	aasdfa	Cua_1	Teo - SeRMN	UsuariNormal
2012-05-24 12:14	2012-05-25 14:55	Normal_7	Cua_1	Teo - SeRMN	UsuariNormal

System managed by **Bumblebee** version 1.3.1, released under the GPL license.
This installation of Bumblebee currently manages 294 users, 57 projects, 17 instruments, 9543 bookings, 2 sample submission forms and 19 requests.
Running under Linux (2.6.32-5-686), Apache (2.2.16), PHP (5.3.3-7+squeeze8, apache2handler mode) with MySQL (5.1.61-0+squeeze1).

Email the **system administrator** for help.

Booking information Copyright © 2012 Servei de Resonància Magnètica Nuclear, UAB

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English ▾

- Main Menu
- Instruments & Queues
- My bookings
- My account
- Change password
- Logout
- IconNMR module
- My samples
- Help contents
- About
- Help

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- **Test** some important information about the queue can be here.
- **Test 2** some important information about the queue can be here.

Sample ID	12
User	296 (UsuariNormal)
Recorded by	296 (UsuariNormal)
Queue	1 (Cua_1)
Sent Date	2012-05-23 18:16
Processed Date	2012-05-25 14:55
My Projects	Teo - SeRMN (T063300 / PK122933 - 2010)

Solvent	AcCapil (Acetone-d6 in Capillary Tube)
Sample name	aasdfa
Amount (mg)	50

Comments

Experiments and properties

Experiment	Teo RMBC
Dangerousness	(Unknown)
Conservation	Temperature between 20-30°C

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Solvent	AcCapil (Acetone-d6 in Capillary Tube)
Sample name	aasdfa
Amount (mg)	50

Comments

Experiments and properties

Experiment (Press CTRL to multiple choice)	[Select one or more options] [No choice] [No choice] [No choice]
Dangerousness (Press CTRL to multiple choice)	[Select one or more options] [No choice] [No choice] Inflammable description
Conservation (Press CTRL to multiple choice)	Temperature between 20-30°C [No choice] [No choice]

Update sample Delete sample Sample as new

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SeRMN Instrument Booking Tool UAB

You are logged as: usuari.normal

- English
- Main Menu
- Instruments & Queues
- My bookings
- My account
- Change password
- Logout
- IconNMR module
- My samples
- Help contents
- About
- Help

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System managed by **Bumblebee** version 1.3.1 (09/03/2012) on a server with 2GB RAM and 10GB disk space. This installation of Bumblebee currently manages 294 users, 57 projects, 17 instruments, 9543 bookings, 2 sample submission forms and 23 requests. Running under Linux (2.6.32-5-686), Apache (2.2.17), MySQL (5.1.61-0+squeezel) and PHP (5.3.17+squeezel, apache2handler mode) with MySQL (5.1.61-0+squeezel1). Email the **system administrator** for help.

Booking information Copyright © 2012 Servei de Resonància Magnètica Nuclear, UAB

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You are logged as: usuari.queue

Si ets gestor d'una cua de rutina...

- English
- Main Menu
- Instruments & Queues
- My bookings
- My account
- Change password
- Logout
- IconNMR module
- My samples
- Manage samples
- Help contents
- About
- Help

Select an instrument to manage your bookings

Instrument	Description	Location
250 MHz - AUTOSERVEI	Bruker Avance DRX-250 autoservei	SeRMN UAB

Select a queue to send a sample - IconNMR

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System managed by **Bumblebee** version 1.3.1 (09/03/2012) on a server with 2GB RAM and 10GB disk space. This installation of Bumblebee currently manages 294 users, 57 projects, 17 instruments, 9543 bookings, 2 sample submission forms and 15 requests. Running under Linux (2.6.32-5-686), Apache (2.2.17), MySQL (5.1.61-0+squeezel) and PHP (5.3.17+squeezel, apache2handler mode) with MySQL (5.1.61-0+squeezel1). Email the **system administrator** for help.

Booking information Copyright © 2012 Servei de Resonància Magnètica Nuclear, UAB

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English ▾

Select an instrument to manage your bookings

- Main Menu
- Instruments & Queues
- My bookings
- My account
- Change password
- Logout
- IconNMR module
- My samples
- Manage samples
- Help contents
- About
- Help

Instrument	Description	Location
250 MHz - AUTOSERVEI	Bruker Avance DRX-250 autoservei	SeRMN UAB

Select a queue to send a sample - IconNMR

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System managed by **Bumblebee** version 1.3.1, released under the GNU GPL. This installation of Bumblebee currently manages 294 users, 57 projects, 17 instruments, 9543 bookings, 2 sample submission forms and 15 requests. Running under Linux (2.6.32-5-686), Apache (2.2.16), PHP (5.3.3-7+squeezed8-apache2-module) with MySQL (5.1.67-0-suse-1.1).

Email the **system administrator** for help.

Booking information Copyright © 2012 Servei de Resonància Magnètica Nuclear, UAB

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Create a iconNMR file

Select the samples you want to s

System managed by **Bumblebee** version 1.3.1, released under the **GNU GPL**.
This installation of Bumblebee currently manages 294 users, 57 projects, 17 instruments, 9543 bookings, 2 sample submission forms and 15 requests.
Running under Linux (2.6.32-5.686), Apache (2.2.16), PHP (5.3.3-7+squeeze8_apache2handler-mod) with MySQL (5.1.61-0+squeeze1)

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My bookings
My account
Change password
Logout

IconNMR module
My samples
Manage samples
Help contents
About
Help

Sent Date	Sample name	User	Queue	Project	Sent by	Select
2012-05-23 11:37	Mostra_Masq_1	usuari.queue	Cua_1	SeRMN - Gestio i Recerca	admin	<input checked="" type="checkbox"/>
2012-05-23 11:38	Queue_2	usuari.queue	Cua_1	Henkel Iberica ATLB	usuari.queue	<input checked="" type="checkbox"/>
2012-05-23 11:38	Queue_1	usuari.queue	Cua_1	Henkel Iberica ATLB	usuari.queue	<input checked="" type="checkbox"/>
2012-05-23 11:40	Normal_1	usuari.normal	Cua_1	Bioiberica	usuari.normal	<input type="checkbox"/>
2012-05-23 11:40	Normal_2	usuari.normal	Cua_1	Bioiberica	usuari.normal	<input checked="" type="checkbox"/>
2012-05-23 11:47	Mostra_2_asnew	usuari.queue	Cua_1	SeRMN - Gestio i Recerca	usuari.queue	<input checked="" type="checkbox"/>
2012-05-23 18:16	aasdfa	usuari.normal	Cua_1	Teo - SeRMN	usuari.normal	<input checked="" type="checkbox"/>
2012-05-24 10:52	sample for nmr	usuari.normal	Cua_1	Teo - SeRMN	usuari.normal	<input type="checkbox"/>
2012-05-24 11:09	another sample for n	usuari.normal	Cua_1	Teo - SeRMN	usuari.normal	<input type="checkbox"/>
2012-05-24 12:14	Normal_7	usuari.normal	Cua_1	Teo - SeRMN	usuari.normal	<input checked="" type="checkbox"/>

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IconNMR module

- My samples
- Manage samples
- Help contents
- About
- Help

File successfully created

Files iconNMR

2012-05-20 14:55 to 2012-05-25 14:55

< earlier | now | later >

Date file	ID file	Queue	Created by
2012-05-25 14:55	2	Cua_1	usuari.queue
2012-05-23 11:58	1	Cua_1	admin

System managed by **Bumblebee** version 1.3.1, released under the **GNU GPL**.
This installation of Bumblebee currently manages 294 users, 57 projects, 17 instruments, 9543 bookings, 2 sample submission forms and 15 requests.
Running under Linux (2.6.32-5-686), Apache (2.2.16), PHP (5.3.3-7+squeeze6, apache2handler mode) with MySQL (5.1.61-0+squeeze1).

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IconNMR module

- My samples
- Manage samples
- Help contents
- About
- Help

Files iconNMR

2012-05-20 14:55 to 2012-05-25 14:55

< earlier | now | later >

Date file	ID file	Queue	Created by
2012-05-25 14:55	2	Cua_1	usuari.queue
2012-05-23 11:58	1	Cua_1	admin

iconNMR files

System managed by **Bumblebee** version 1.3.1, released under the **GNU GPL**.
This installation of Bumblebee currently manages 294 users, 57 projects, 17 instruments, 9543 bookings, 2 sample submission forms and 15 requests.
Running under Linux (2.6.32-5-686), Apache (2.2.16), PHP (5.3.3-7+squeeze8, apache2handler mode) with MySQL (5.1.61-0+squeeze1).

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What should Iceweasel do with this file?

Open with gedit (default)

Save File

Do this automatically for files like this from now on.

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System managed by **Bumblebee** version 1.3.1, released under the GPL v2.
 This installation of Bumblebee currently manages 294 users, 57 projects, 17 instruments, 9543 bookings, 2 sample submission forms and 15 requests.
 Running under Linux (2.6.32-5-686), Apache (2.2.16), PHP (5.3.3-7+squeeze8, apache2handler mode) with MySQL (5.1.61-0+squeeze1).

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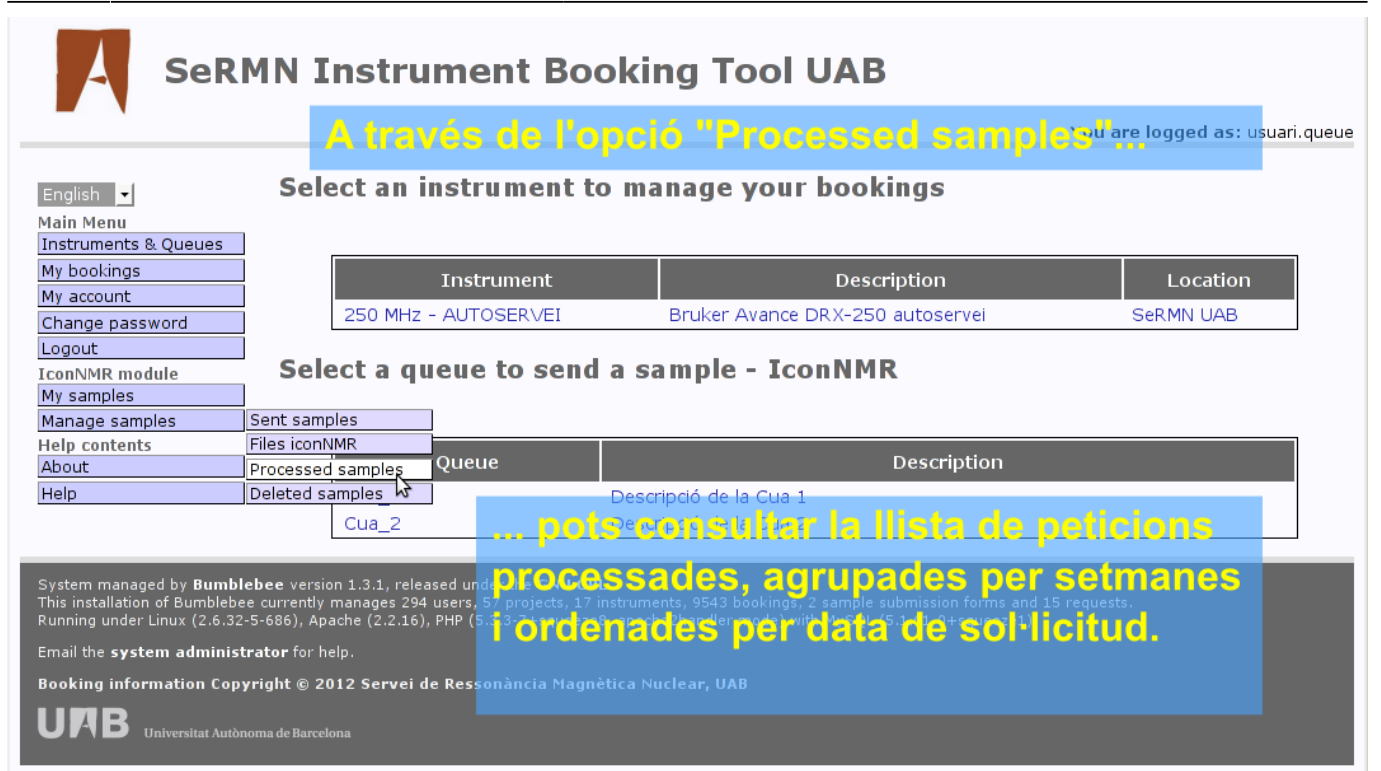
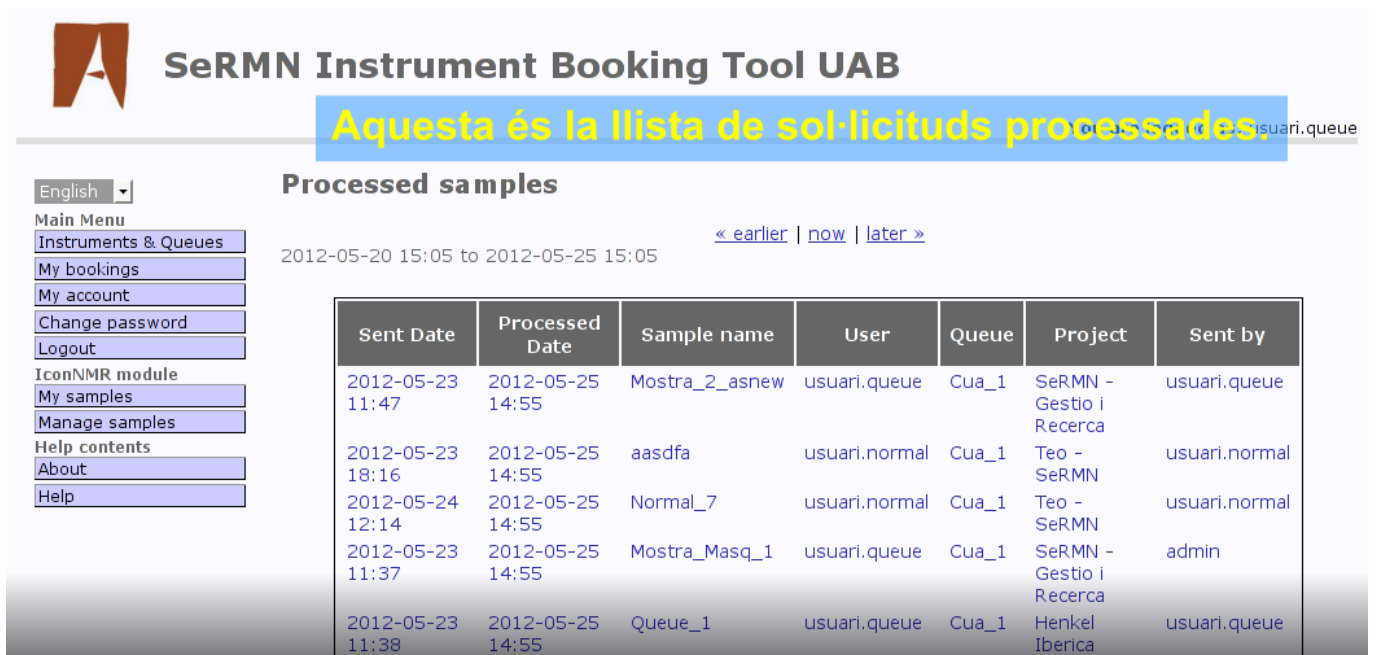
System managed by **Bumblebee** version 1.3.1, released under the GPL license.
This installation of Bumblebee currently manages 294 users, 57 projects, 17 instruments, 9543 bookings, 2 sample submission forms and 15 requests.
Running under Linux (2.6.32-5-686), Apache (2.2.16), PHP (5.3.3-7) and the Bump database backend mode with MySQL (5.1.41+gucos3).

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English

Main Menu

- Instruments & Queues
- My bookings
- My account
- Change password
- Logout
- IconNMR module
- My samples
- Manage samples
- Help contents
- About
- Help

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Sample ID	12
User	296 (UsuariNormal)
Recorded by	296 (UsuariNormal)
Queue	1 (Cua_1)
Sent Date	2012-05-23 18:16
Processed Date	2012-05-25 14:55
All Projects	Teo - SeRMN (T063300 / PK122933 - 2010-10-01)

Solvent	AcCapil (Acetone-d6 in Capillary Tubes)
Sample name	aasdfa
Amount (mg)	50

Comments

Experiments and properties

Experiment	Teo HMBC Teo Proto
Dangerousness	(Unknown)
Conservation	Temperature between 20-30°C

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